
HB-1D

Hybridiser
Operator's Manual

Issue 21.
Date of issue:
4th March 2003
® The Techne logo is the trademark
of Techne and Techne Inc.
© Techne, 2003

CONTENTS

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HB-1D issue 20 onwards Operator's Manual	page
SAFETY AND INSTALLATION	
Declaration of Conformity	3
English	4
Français	5
Deutsch	6
Español	7
Italia	8
INTRODUCTION	9
Initial checks	9
Specification	10
Unpacking	11
Installation	11
Warning	13
Servicing by Techne and their Agents	13
Switching on	14
Temperature control	14
Speed control	14
OPERATION	
The tubes	15
The HB-1D	17
The front panel controls	18
Basic Protocol for Southern Hybridisations	20
When the Protocol is Finished	20
TECHNICAL INFORMATION	21
Operator Maintenance	21
Accessories and Spares	23

Techne
Duxford
Cambridge
CB2 4PZ

Declaration of Conformity

Techne's HB-1D complies with the following European Standard:

EN 61326:1998 Electrical equipment for measurement, control and laboratory use.
EMC requirements.

It has also been designed to comply with the following European Standards:

EN 61010-1:1993 Safety requirements for electrical equipment for measurement, control and laboratory use.

EN 61010-2-010:1995 Particular requirements for laboratory equipment for the heating of materials.

I have made all reasonable enquiries regarding the unit stated and its conformance to the following EU directives:

Low Voltage directive, 73/23/EEC and amendment 93/68/EEC, and

EMC Directive 89/336/EEC and amendments 91/263/EEC and 92/31/EEC.

To the best of my knowledge and belief these units conform to these directives.

Techne's HB-1D also complies with CFR 47 Part 15B for a class A digital device.

CFR 47:1997 Code of Federal Regulations: Part 15 Subpart B-Radio Frequency Devices - Unintentional Radiators

This Declaration is controlled under an ISO 9001:1994 system certificated by BSI Quality Assurance, certificate number FM13585.

Signature

Name

B C Coombes

Position

Quality Manager

Issue 20

18/11/2002

Introduction

Please read all the information in this booklet before using the unit.

Warning

HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material.

Techne have taken great care in the design of these units to protect operators from hazards, but users should pay attention to the following points:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS;
- DO NOT put hot objects on or near combustible objects;
- DO NOT operate the unit close to inflammable liquids or gases;
- DO NOT place any liquid directly in your unit;
- At all times USE COMMON SENSE.

Operator Safety

All users of Techne equipment must have available the relevant literature needed to ensure their safety.

It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Techne the protection provided by the equipment to the user may be impaired.

All Techne units have been designed to conform to international safety requirements and are fitted with an overtemperature cutout. The cutout is preset to protect the unit.

If a safety problem should be encountered, switch off at the mains socket and remove the plug from the supply.

Installation

1. All Techne units are supplied with a power cable. This may be integral or plug-in.
2. Before connecting the mains supply, check the voltage against the rating plate. Connect the mains cable to a suitable plug according to the table below. **Note that the unit must be earthed to ensure proper electrical safety.**

Connections	220/240V	110/120V
Live	Brown	Black
Neutral	Blue	White
Earth	Green/yellow	Green

The fused plug supplied with the mains lead for use in the UK is fitted with the following value fuse to protect the cable:

230V UK	5 AMP
---------	-------

The fuse in the unit protects the unit and the operator.

Note that units marked 230V on the rating plate work at 220V; units marked 120V work at 110V. In both cases, however, the heating rate will degrade by approximately 8%. The rating plate is on the rear of the unit.

3. Plug the mains cable into the socket on the rear of the unit.
4. Place the unit on a suitable bench or flat workspace, or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction.
5. Symbols on or near the power switch of the unit have the following meanings:

I	: mains switch On
O	: mains switch Off

After use

When you have finished heating samples, remember that parts of the unit – the tubes, blocks and associated accessories – may be very hot. Take the precautions listed earlier.

Guarantee

The unit is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is from the date of purchase, and within this period all defective parts will be replaced free of charge provided that the defect is not the result of misuse, accident or negligence. Servicing under this guarantee should be obtained from the supplier.

Notwithstanding the description and specification(s) of the units contained in the Operator's Manual, Techne hereby reserves the right to make such changes as it sees fit to the units or to any component of the units.

This Manual has been prepared solely for the convenience of Techne customers and nothing in this Instruction Book shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the units or components.

User maintenance

NOTE THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL. REMOVING THE INSIDE FRONT OR REAR PANELS EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES. THERE ARE NO USER MAINTAINABLE PARTS WITHIN THE EQUIPMENT.

In the unlikely event that you experience any problems with your unit which cannot easily be remedied, you should contact your supplier and return the unit if necessary. Please include any details of the fault observed and remember to return the unit in its original packing. Techne accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier. See the Decontamination Certificate supplied with your unit.

1. Cleaning

Before cleaning your unit ALWAYS disconnect from the power supply and allow to cool below 50° C.

Your unit can be cleaned by wiping with a damp soapy cloth. Care should be exercised to prevent water from running inside the unit. Do not use abrasive cleaners.

2. Overtemperature cutout

The overtemperature cutout may be a sensitive mechanical device and mechanical shock can cause it to trip.

- In the event of no heater power, check the mains plug and lead, then reset the cutout control.
- Repeated operation of the cutout indicates a serious fault: you may need to return the unit to your supplier for repair.

3. Fuses

Your unit is protected by one or two fuses. These should only be changed by suitably qualified personnel.

If the fuses blow persistently, a serious fault is indicated and you may need to return the unit to your supplier for repair.

Contact Information

For technical, sales or servicing information, contact your local Techne dealer or,

Techne, Duxford,

CAMBRIDGE, CB2 4PZ, United Kingdom.

Telephone: 01223 832401

Telefax: 01223 836838

Service: 01223 836950

Out of office hours, telephone or fax

e-mail: sales@techne.com

Web site: www.techne.com

or,

Techne Incorporated,

University Park Plaza, 743 Alexander Road,

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INTRODUCTION

Read the whole of this book before commencing work with the unit.

Blotting techniques in which DNA, RNA or protein are immobilised onto nylon or nitrocellulose filters form a cornerstone of molecular biology. Detection of the desired sequence is usually performed using radioactive probes. Given the expense and hazardous nature of the probes, probing should:

- a) be performed under safe conditions
- b) use minimal volumes (to improve signal and reduce expense)
- c) allow reuse of materials
- d) be simple to use

The Techne HB-1D meets all these requirements.

Up to five 25 x 25cm filters can be probed in one of the large hybridisation tubes. Hybridisations can be performed in a minimal volume of 5ml, an 80% reduction over conventional systems, reducing expense and improving signal. The operator is protected from irradiation throughout and the apparatus is extremely simple to use. However, the chemical manufacturer's safety precautions must be used when using the HB-1D.

The Techne HB-1D also incorporates the following features:

- a) a temperature range of ambient +10°C to 100°C.
- b) it holds 1, 2, 3 or 4 x 80mm diameter hybridisation tubes.
- c) it will hold 3, 6, 9 or 12 x 44mm diameter tubes if you use a three-tube carrier.
- d) it will hold 4, 8, 12 or 16 mini tubes if you use a four-tube carrier.
- e) the HB-1D comes with adjustable feet to enable accurate levelling.

Initial Checks

When you receive your new HB-1D remove all the packing. Check the unit for any transit damage. If there is any transit damage at all inform Techne Sales Department or your Techne dealer immediately.

When unpacking the unit, check that the following have been removed from the packing: **a power cable; a tube of silicon grease; a decontamination certificate and a Guarantee Card.**

Check that the unit is the correct rating for your power supply by looking on the Rating Plate which is on the rear of the unit.

Within the guarantee period, shown on the Guarantee Card, we undertake to supply replacements free of charge for parts which may, on examination, prove to be defective, provided that the defect is not the result of misuse, accident or negligence. Any instrument requiring service under this guarantee should be sent to the supplier through whom it was purchased, or, in the case of difficulty, it should be carefully packed in its original packing and consigned, carriage paid, to us. Techne takes no responsibility for returned goods damaged in transit.

Returned goods will not be processed without a Returns Authorisation Number. Call + (44) (0)1223 832401 for a number. On all correspondence, please quote the Serial Number in full and/or the Sales Order Number. **Please write the Returns Number on the outside of any packing.**

Specification

Maximum Number of Hybridisation Tubes	6 (large and small) 24 mini tubes
Temperature Range #	ambient +10° to 100°C (NOTE: minimum temperature must be greater than 10°C above ambient)
Rotation of Hybridisation Tubes	5 rpm to 20rpm and off
4 steps, 5 positions and off	Approx steps of 4 rpm
Adjustable rocking platform	5-20 or 15-60 opm
Temperature stability in chamber	better than 1.0°C
Stability in the tubes	better than +0.1°C
Temperature uniformity in the chamber *	better than +1.5°C
Temperature uniformity in the tubes *	better than +1.0°C
Temperature set point resolution	0.1°C
Absolute accuracy	better than +0.3°C
Warm Up Time Within Tube (16° to 100°C)	50 mins
Heater Power	750 W
Temperature selection	Digital
Temperature Display	LED
Temperature Sensor	PRT
External Dimensions	
Width	385mm
Depth	285mm
Height	555mm
Internal Dimensions	
Width	260mm
Depth	210mm
Height	330mm
Weight	
Net	20Kg

The Hybridiser includes a preset re-settable overtemperature cutout.

* Temperature measured in the centre of each tube.

Large tubes are fitted with 70°C seals, seals part number 6101488 are required to work to 100°C.

For the Hybridisation Tubes see the list under Accessories

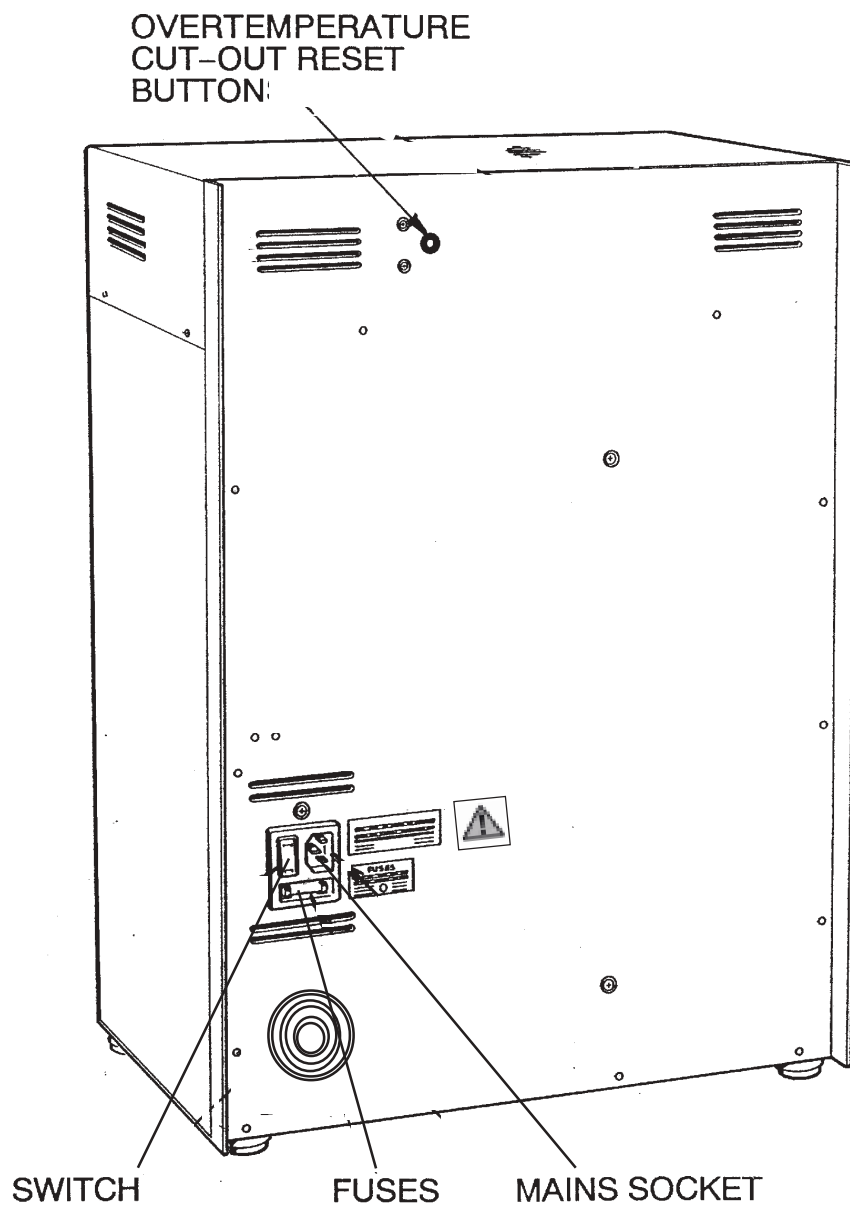
Unpacking

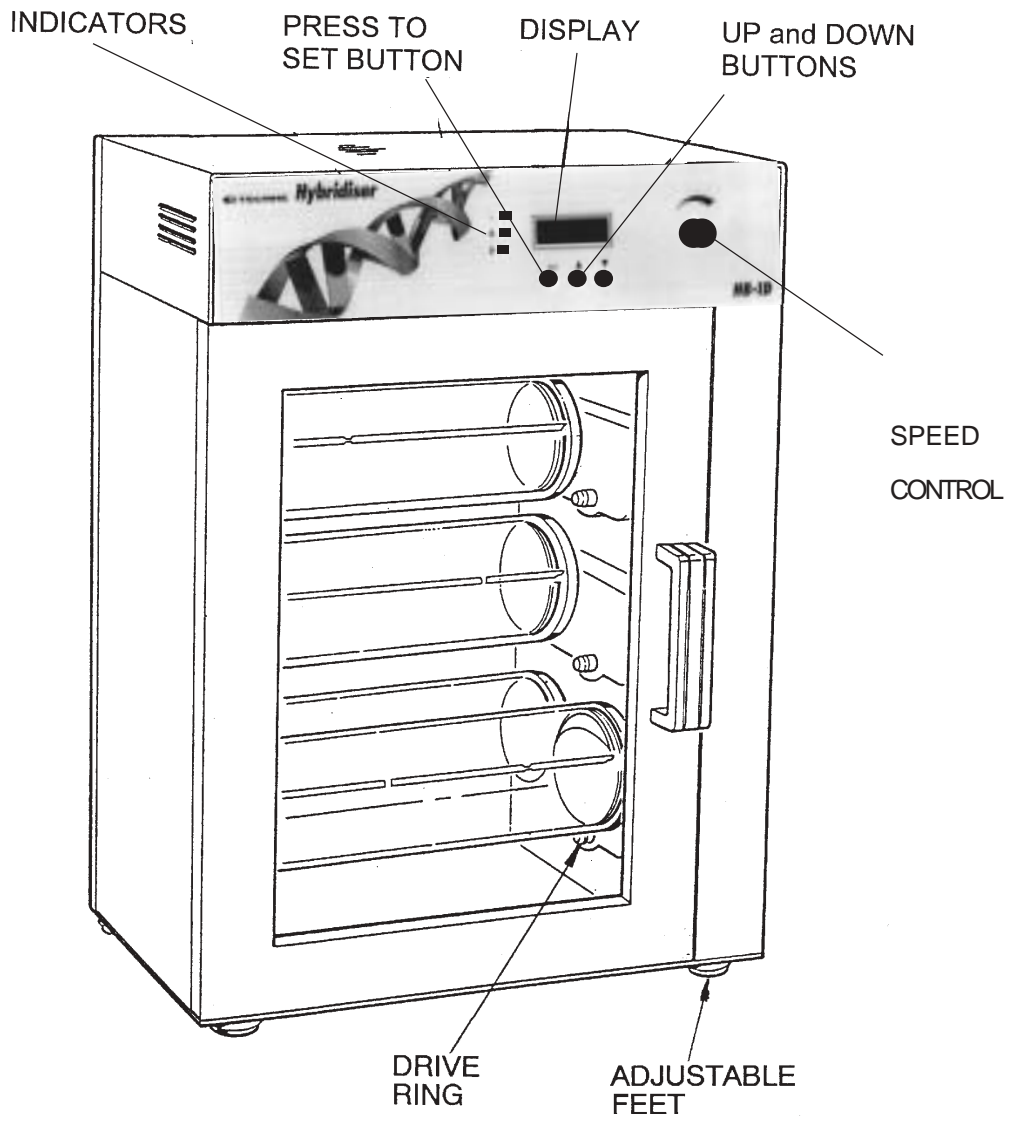
Ensure that the following have been removed from the packing:

- The Hybridiser.
- Drip Tray (not Techne Inc).
- Power Lead.
- Guarantee Card.
- Decontamination Certificate.
- 3 Hybridisation tubes complete with end caps, seals and grease.
- Extra Tubes if ordered especially.

Installation

- 1 Place the unit on a suitable bench, a flat workspace, or in a safety cupboard if required, ensuring that the air inlet and outlet louvres are not obstructed.
- 2 Insert some liquid into a tube and place the tube in the Hybridiser. Adjust the feet by unscrewing them anticlockwise slightly until the unit is level. Check the level by putting the tube in several positions.
- 3 Plug the mains cable into the socket on the rear of the unit.

Warning

THE UNIT WORKS UP TO 100°C. This temperature can cause burns. Techne take great efforts to protect operators from hazards, but the following points are brought to your attention:

- 1 Use care and wear protective gloves to remove hybridisation tubes from the cabinet if the unit is at high temperature.

If the protocol allows, it is better to allow the unit to cool to below 40°C before attempting to remove the tubes.
- 2 The coloured end cap of the 80mm large tube and the 44mm small tube has two holes, so that you have to be careful not to spill the probe through these.

One is the hole which allows the air to expand when the tube is heated.

The other is the filling hole which is revealed when the spindle cap is removed.

DO NOT TIP THE TUBE SO THAT THE COLOURED END IS DOWN AS THIS WILL CAUSE A SPILLAGE.
- 3 The unit has been designed to conform to international safety specifications and is fitted with a resettable over-temperature cutout and an adjustable cutout. If, for any reason, the temperature rises above 115°C, the over-temperature protection cuts out the heater, but the tubes continue to rotate. The over-temperature light comes on. Check the reason for the cut out and, if there is no obvious cause, reset the cutouts once the unit has cooled. The reset, on the back of the unit, is operated by pressing the centre of the button indicated on page 11.

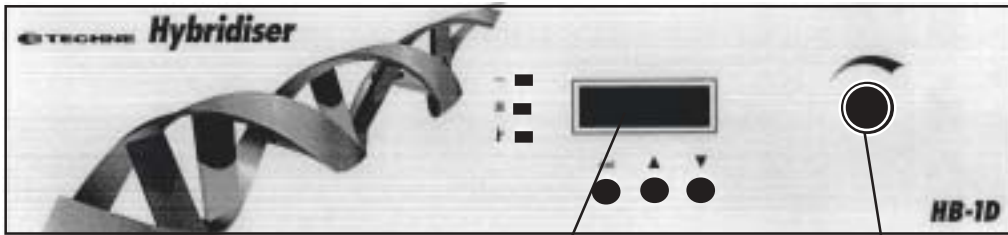
Servicing by Techne and their Agents

Please note that neither Techne nor any of their agents can accept the return of ANY GLASSWARE from these units.

If the unit is returned it must be thoroughly decontaminated and a CERTIFICATE OF CLEANLINESS returned with the unit.

Switching On

Once your HB-1D is connected to a power supply, switch it on using the power switch next to the power inlet socket on the rear panel.


Temperature display

Speed control knob

Front panel of the HB-1D

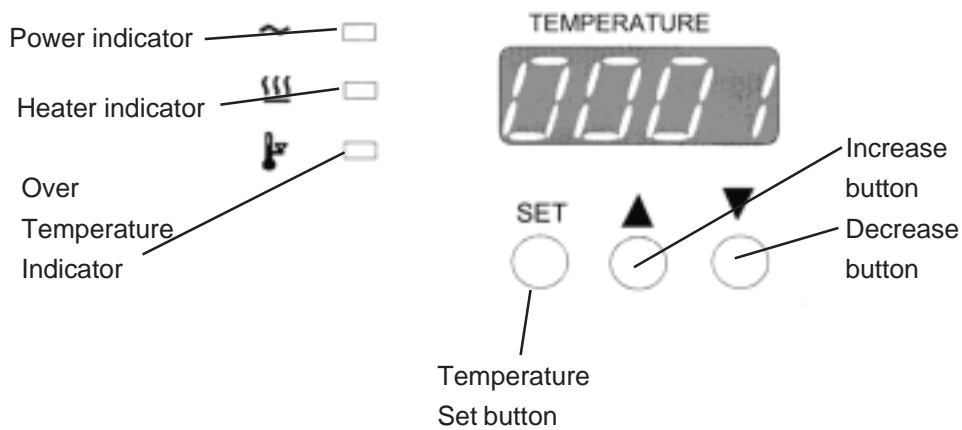Temperature Control

Controller Display

When you first switch on, the display will show the edition of the software which your unit uses. For example software issue "1" would be shown as follows:

It will display this for 1 second, then the actual temperature of the incubator will be indicated.

Speed control

Fully anticlockwise the speed control knob switches the rotation of the tubes off

It then increases the speed clockwise and decreases the speed anticlockwise in steps of approximately 4 rpm. Switch the knob until you have the speed you require.

OPERATION

The Tubes

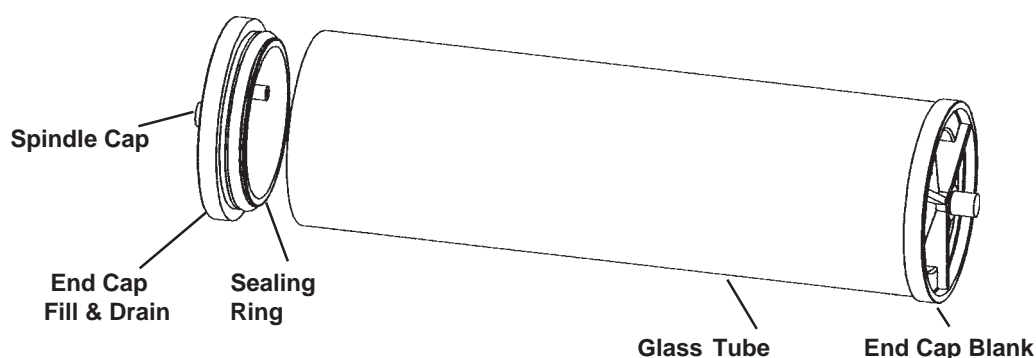
The Hybridisation Tubes must be checked regularly and replaced if any damage occurs. Any damage to the ends of the tubes can cause them to break when the end caps are inserted.

Each tube must be fitted with a radiation shielding end cap if you are using a radioactive probe. These end caps are supplied with all tubes which are supplied with the HB-1D and any other tubes supplied after December 1999. Normal handling procedures must be observed when handling any tube.

Load the membrane and reagents according to which size hybridisation tube you are using.


Large tube (80mm diameter) assembly

Remove the tube end cap which has the holes in it. Place the filters in the tube and replace the end cap. A small quantity of Techne O-ring grease, see the list of accessories, **must** be smeared on the seal to facilitate the insertion of the end cap. Remove the spindle cap and the liquids appropriate to the protocol may then be added through the hole in the spindle using a syringe. Replace the spindle cap and press it onto the spindle so that it clicks into place. Note: the tube will not rotate properly if the cap is not fitted securely

Mini tube (4 x 32mm dia) assembly

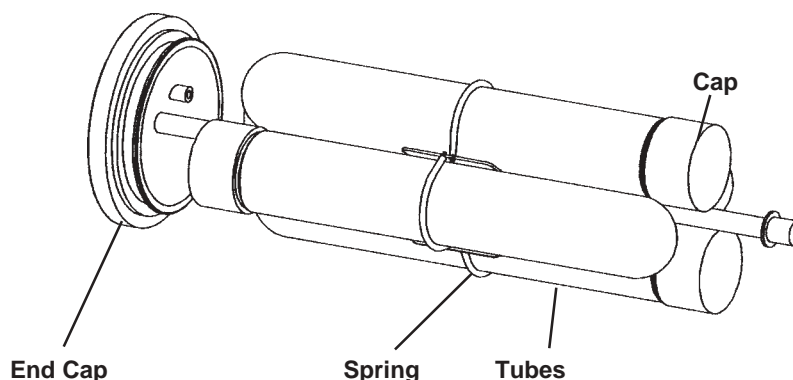
Unscrew the end cap of the tube. Place the filters and the liquids appropriate to the protocol in the tube and replace the end cap. Fix the tube into a multi-tube holder, using the springs, and ensure that there are four tubes in the holder. The end caps must be as shown, on alternate ends. The multi-tube holder will not rotate properly if there are not four mini tubes in it.

Small Tube (3 x 44mm dia) assembly

Unscrew the tube end cap. Place the filters in the tube and replace the end cap. Remove the spindle cap and the liquids appropriate to the protocol may then be added through the hole in the spindle using a syringe. Replace the spindle cap and press it onto the spindle so that it clicks into place.

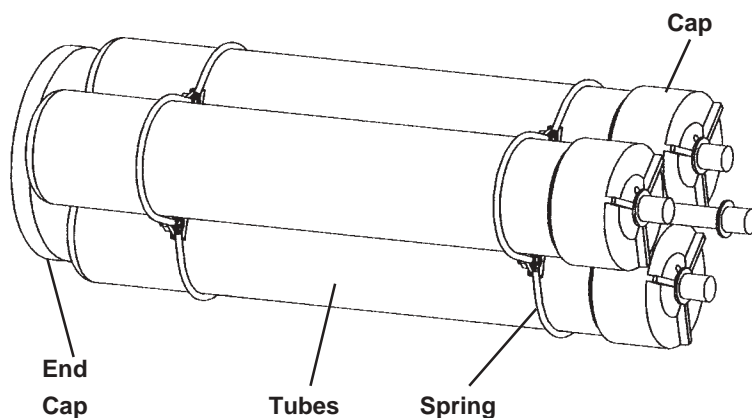
The assembly is similar to the four tube assembly but the end caps do not have to be at different ends of the carrier. Fix the tube into a multi-tube holder, using the springs, and ensure that there are three tubes in the holder. The multi-tube holder will not rotate properly if there are not three small tubes in it.

16 x 15ml Tube assembly

Unscrew the end cap from the tube. Place the filters and the liquids appropriate to the protocol in the tube and replace the end cap.

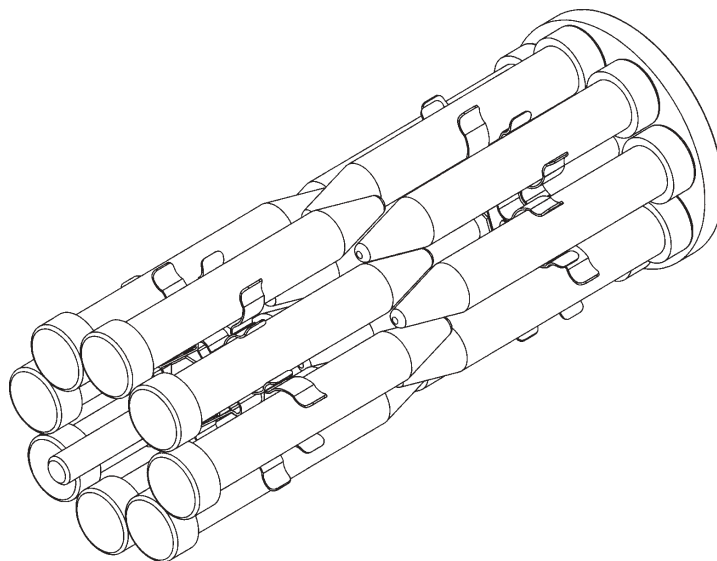Assemble the tubes into a multi-tube holder, using the spring clips. The multi-tube holder will not rotate properly if it is not balanced properly. Preferably the multi-tube holder should be filled completely even if all the tubes are not being used.


The HB-1D

- 1 Insert the tubes or tube holder into the HB-1D ensuring that they rotate freely and engage on the drive rollers.
- 2 Close the door.
- 3 Connect the mains electricity supply to the mains socket at the rear of the unit. Switch the power ON from the switch located next to the mains input lead. The POWER indicator and the display on the front panel will light.
- 4 The Heater Indicator will be lit while the heater is on.

The Front Panel Controls


The front panel controls consist of three buttons for controlling the display, a four digit LED display and three indicators.

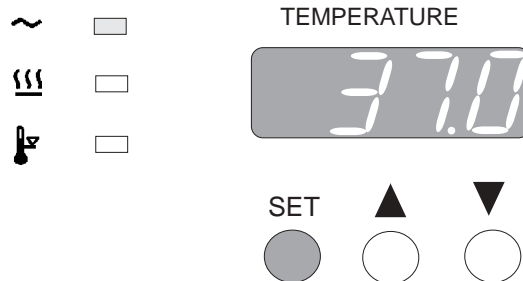The SET temperature Button


The **SET** temperature button displays the set temperature when pressed.

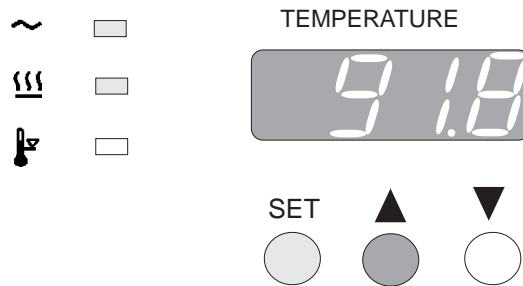The UP ARROW Button

When the **SET** temperature button is held down and the **UP ARROW** button is pressed, the set temperature is increased.

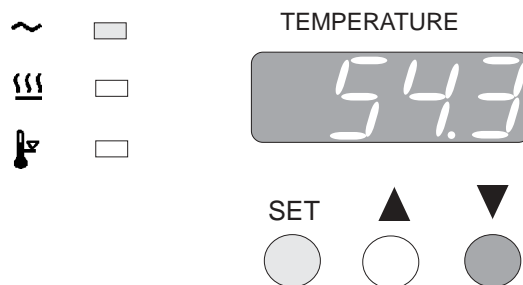The DOWN ARROW Button

When the **SET** temperature button is held down and the **DOWN ARROW** button is pressed, the set temperature is decreased.

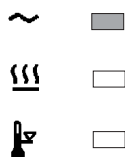
Speed of Change of Set Temperature

Each press of the **UP ARROW** or **DOWN ARROW** buttons will increase or decrease the set temperature by .1°C. If the buttons are held down the temperature change will accelerate to 5° per second.

Tube temperature

NOTE: It takes approximately 50 minutes for the interior of the hybridisation tubes to reach the displayed temperature depending upon the number of filters and amount of hybridisation fluid used.

Power Indicator

TEMPERATURE

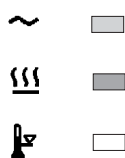
SET

The top indicator shows that there is power to the unit

Heater Indicator

TEMPERATURE

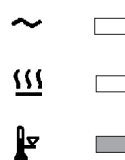
SET

The next indicator shows when the heater is heating the unit. When the temperature is being set, and the new set temperature is higher than the temperature already in the unit, the heater indicator will light as the unit tries to follow the set temperature. If the light is on continuously the heater is getting constant power. The only exception is described under Over-Temperature Indicator. As the temperature approaches the set temperature the heater indicator will flash. When set temperature is reached the indicator will stay on for shorter periods. If the oven is above the set temperature then the indicator will be off, as the heater is not getting any power.

Over-Temperature Indicator

TEMPERATURE

SET

If the unit should, for any reason, exceed the temperature set for the over-temperature cutout, the over-temperature indicator will light. The heater will have been switched off and the unit will begin to cool even if the heater light is on (the light staying on or not depends on which circuit has sensed an over-temperature). The fan will continue to circulate the air.

BASIC PROTOCOL FOR SOUTHERN HYBRIDISATIONS *

INSERTION OF FILTERS

Place a wet or moist filter, DNA side facing inwards, into a hybridisation tube and allow it to adhere to the glass. Large trapped air bubbles should be smoothed out with a glass pipette; smaller air bubbles are acceptable and do not interfere with hybridisation so long as the surface of the filter is covered with hybridisation fluid.

Several filters may be placed in one tube; there is a slight loss of signal, but no increase in background in the lower filters. To improve the signal we recommend the use of separation membranes, part number FMEM2020.

PRE-HYBRIDISATION

Add 5 or 10ml of fluid, according to the number of filters and the size of tube, so that the filters are covered. Allow to pre-hybridise for usual period.

HYBRIDISATION

Drain off the pre-hybridisation fluid by inverting the tube. Add hybridisation liquid with radioactive probe added by using a syringe and needle through the hole in the blue end-cap. The minimum volume for a single filter is 5ml.

WASHES

Washing depends on the stringency required.

One method is to wash genomic blots in the following fashion:

- 1) 2 x SSC, 50ml for 10 minutes
- 2) 2 x SSC, 50ml for 10 minutes
- 3) 0.1 x SSC, 0.1% SDS x 50ml for 10 minutes

This gives very clean backgrounds.

NOTE

- 1) All radioactive waste can be collected for disposal; none need be disposed of down the sink.
- 2) Radioactive probes can be collected and reused several times.
- 3) As the protocols for this unit usually involve the use of radioactive materials the appropriate manufacturer's safety precautions must be taken into consideration.

* Courtesy of: Dr M J S Dyer.

When the Protocol is Finished

When the protocol has finished switch off the unit. Open the door and remove the tubes, again being careful to keep the coloured end-cap uppermost. Remember the tubes may be hot. Either wear gloves or, if the protocol permits, allow the tubes to cool before removal.

TECHNICAL INFORMATION

Brief fault finding notes are given in this section.

NOTE THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL. REMOVING THE CASE EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES. THERE ARE NO OPERATOR MAINTAINABLE PARTS WITHIN THIS EQUIPMENT.

Operator Maintenance

General

Should you experience any problems with your HB-1D which cannot be easily remedied, you should contact your supplier and return the unit if necessary.

Fill in the Decontamination Certificate provided with the unit and return it to Techne Service department giving as much information as possible to help diagnose the fault. Please remember to return the unit in its original packing, Techne accept no responsibility for damage to units which are not properly packed: if in doubt, contact your supplier.

Overtemperature Cutout

The overtemperature cutout is a sensitive mechanical device and mechanical shock can cause it to trip.

In the event of no heater power, check the mains plug and lead, then reset the cutout control by depressing the centre of the button.

Repeated operation of the cutout indicates a serious fault: you may need to return your unit to your supplier for repair.

Fuses

If the power indicator fails to light when power is supplied to the unit the fuses may have blown. Check that there is no external cause such as a faulty plug; check both fuses and replace any faulty fuse with one of the correct value. Fuse values are given on the label next to the mains input socket.

O ring seals on the Hybridisation Tubes

The O ring seals on the 80mm dia Hybridisation Tubes must be checked regularly and replaced before a leak occurs. Each tube is provided with 6 spare O rings and further O rings can be obtained from Techne or your dealer.

Hybridisation Tubes

The Hybridisation Tubes must be checked regularly and replaced if any damage occurs. Any damage to the ends of the tubes can cause them to break when the end caps are inserted. Further Tubes can be obtained from Techne or your dealer.

Cleaning your HB-1D

Before cleaning your unit ALWAYS disconnect from the power supply and allow to cool.

You can clean the case of the HB-1D with a cloth dampened in soapy water or ethanol. No part of the unit should be immersed in solvents. Do not use acetone or abrasive cleaners.

If You Spill the Fluid

Because of the nature of operation of the HB-1D, the complete ductwork system must be cleaned if a spillage does occur. This is to remove any airborne particles which may have been blown around the system and to guarantee full operator safety.

Switch the HB-1D off and disconnect it from the power supply. Follow this sequence of operations when the HB-1D has been allowed to cool below at least 50°C:

- a) Remove all glassware from the HB-1D.
- b) Remove the drip tray.
- c) Clean all surfaces of the inner chamber using a suitable cleaning and decontaminating solution.

Be careful of the drive shafts.

Cleaning and decontamination materials which Techne recommend include:

Decon 90

Neutracon

or mild detergent.

If an Incorrect Message is Displayed

If a message is displayed which shows a strange temperature or HELP, the PRT (temperature sensor) has probably gone open circuit. This could mean that a wire has broken; a wire or the connector to the PCB has become disconnected; or the PRT has been broken somehow; etc. You should contact your supplier and return the unit if necessary. Remember to fill in the Decontamination Certificate provided with the unit and return it to Techne Service department or your dealer giving as much information as possible to help diagnose the problem.

Accessories and Spares

The following may be purchased from Techne or their dealers. All Techne tubes are made from Borosilicate glass. These tubes are the preferred tubes:

FHB12	80mm external diameter hybridisation tube assembly complete with sealing rings and end caps
FHB11	80mm external diameter tube only
6101082	End cap with holes (blue) less sealing ring
6101387	Spindle cap (for use with 6101082).
6101083	End cap without holes (natural) less sealing ring
6101090	Sealing ring (for 70°C operation)
6101488	Sealing ring (accessory for 100°C operation)
6101351	Techne O-ring Grease (tube)
FHB32	44mm external diameter hybridisation tube assembly complete with seal and screw on end cap
FHB31	44mm external diameter tube only
6007625	Seal (for use with 44mm external diameter tubes).
6101486	Screw on end cap less seal.
6101387	Spindle cap (for use with 6101486).
FHB42	4 x 32mm external diameter hybridisation tubes complete with screw caps, seals and multi-tube holder
FHB41	4 x 31mm external diameter tubes only
FHB43	Multi-tube holder
6102375	End cap for the mini tube
6102376	Seal for the mini tube
6101387	Spindle cap
6101828	Spring for the multi-tube holder
FHBSH1	Shelf
FHBST1	Rocking platform 20/40 Oscillations
FHBST2	Rocking Platform 20/60 Oscillations
FHTRACK	Hybridisation Tube Rack
FMEM2020	Separation Membrane 20cm x 20cm

Contact Techne Sales or your dealer for further details.