Hybrigene

Hybridiser
Operator's Manual

Issue 2.
Date of issue:
27th February 2003
® The Techne logo is the trademark
of Techne and Techne Inc.
© Techne, 2003

CONTENTS

Hybrigene Operator's Manual	page
SAFETY AND INSTALLATION	
Declaration of Conformity	3
English	4
Français	5
Deutsch	6
Español	7
Italia	8
INTRODUCTION	9
Initial checks	9
Specification	10
Installation	12
Fitting stacking plates and wall plates	13
Fixing procedures	14
Switching on	15
Temperature control	15
Warning	16
Servicing by Techne and their Agents	16
OPERATION	
The tubes	17
The Hybrigene	19
Basic Protocol for Southern Hybridisations	20
When the Protocol is Finished	21
TECHNICAL INFORMATION	22
Operator Maintenance	22
Accessories and Spares	23

Techne
Duxford
Cambridge
CB2 4PZ

Declaration of Conformity

Techne's Hybrigene complies with the following European Standard:

EN 61326:1998 Electrical equipment for measurement, control and laboratory use.
EMC requirements.

It has also been designed to comply with the following European Standards:

EN 61010-1:1993 Safety requirements for electrical equipment for measurement, control and laboratory use.

EN 61010-2-010:1995 Particular requirements for laboratory equipment for the heating of materials.

I have made all reasonable enquiries regarding the unit stated and its conformance to the following EU directives:

Low Voltage directive, 73/23/EEC and amendment 93/68/EEC, and

EMC Directive 89/336/EEC and amendments 91/263/EEC and 92/31/EEC.

To the best of my knowledge and belief these units conform to these directives.

Techne's Hybrigene also complies with CFR 47 Part 15B for a class A digital device.

CFR 47:1997 Code of Federal Regulations: Part 15 Subpart B-Radio Frequency Devices - Unintentional Radiators

This Declaration is controlled under an ISO 9001:1994 system certificated by BSI Quality Assurance, certificate number FM13585.

Signature	
Name	B C Coombes
Position	Quality Manager
Issue 1	01/08/2002

Introduction

Please read all the information in this booklet before using the unit.

Warning

HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material.

Techne have taken great care in the design of these units to protect operators from hazards, but users should pay attention to the following points:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS;
- DO NOT put hot objects on or near combustible objects;
- DO NOT operate the unit close to inflammable liquids or gases;
- DO NOT place any liquid directly in your unit;
- At all times USE COMMON SENSE.

Operator Safety

All users of Techne equipment must have available the relevant literature needed to ensure their safety.

It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Techne the protection provided by the equipment to the user may be impaired.

All Techne units have been designed to conform to international safety requirements and are fitted with an overtemperature cutout. The cutout is preset to protect the unit.

If a safety problem should be encountered, switch off at the mains socket and remove the plug from the supply.

Installation

1. All Techne units are supplied with a power cable. This may be integral or plug-in.
2. Before connecting the mains supply, check the voltage against the rating plate. Connect the mains cable to a suitable plug according to the table below. **Note that the unit must be earthed to ensure proper electrical safety.**

<i>Connections</i>	<i>220/240V</i>	<i>110/120V</i>
Live	Brown	Black
Neutral	Blue	White
Earth	Green/yellow	Green

The fused plug supplied with the mains lead for use in the UK is fitted with the following value fuse to protect the cable:

230V UK 5 AMP

The fuse in the unit protects the unit and the operator.

Note that units marked 230V on the rating plate work at 220V; units marked 120V work at 110V. In both cases, however, the heating rate will degrade by approximately 8%. The rating plate is on the rear of the unit.

3. Plug the mains cable into the socket on the rear of the unit.
4. Place the unit on a suitable bench or flat workspace, or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction.
5. Symbols on or near the power switch of the unit have the following meanings:

I : mains switch On
 O : mains switch Off

After use

When you have finished heating samples, remember that parts of the unit – the tubes, blocks and associated accessories – may be very hot. Take the precautions listed earlier.

Guarantee

The unit is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is from the date of purchase, and within this period all defective parts will be replaced free of charge provided that the defect is not the result of misuse, accident or negligence. Servicing under this guarantee should be obtained from the supplier.

Notwithstanding the description and specification(s) of the units contained in the Operator's Manual, Techne hereby reserves the right to make such changes as it sees fit to the units or to any component of the units.

This Manual has been prepared solely for the convenience of Techne customers and nothing in this Instruction Book shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the units or components.

User maintenance

NOTE THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL. REMOVING THE INSIDE FRONT OR REAR PANELS EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES. THERE ARE NO USER MAINTAINABLE PARTS WITHIN THE EQUIPMENT.

In the unlikely event that you experience any problems with your unit which cannot easily be remedied, you should contact your supplier and return the unit if necessary. Please include any details of the fault observed and remember to return the unit in its original packing. Techne accept no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier. See the Decontamination Certificate supplied with your unit.

1. Cleaning

Before cleaning your unit ALWAYS disconnect from the power supply and allow to cool below 50° C. Your unit can be cleaned by wiping with a damp soapy cloth. Care should be exercised to prevent water from running inside the unit. Do not use abrasive cleaners.

2. Overtemperature cutout

The overtemperature cutout may be a sensitive mechanical device and mechanical shock can cause it to trip.

- In the event of no heater power, check the mains plug and lead, then reset the cutout control.
- Repeated operation of the cutout indicates a serious fault: you may need to return the unit to your supplier for repair.

3. Fuses

Your unit is protected by one or two fuses. These should only be changed by suitably qualified personnel. If the fuses blow persistently, a serious fault is indicated and you may need to return the unit to your supplier for repair.

Contact Information

For technical, sales or servicing information, contact your local Techne dealer or,

Techne, Duxford,
 CAMBRIDGE, CB2 4PZ, United Kingdom.
 Telephone: 01223 832401
 Telefax: 01223 836838
 Service: 01223 836950
 Out of office hours, telephone or fax
 e-mail: sales@techne.com
 Web site: www.techne.com

or,
 Techne Incorporated,
 University Park Plaza, 743 Alexander Road,
 Princeton, New Jersey, 08540-6328, USA.
 Telephone: (609) 452-9275
 Toll free: 1-800-225-9243
 Telefax: (609) 987-8177
 e-mail: salesus@techne.com
 Web site: www.techneusa.com

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Techne, Duxford,
CAMBRIDGE, CB2 4PZ, United Kingdom.

Telefono: +44 (0)1223 832401

Telefax: +44 (0)1223 836838

Assistenza: +44 (0)1223 836950

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e-mail: sales@techne.com

Web site: www.techne.com

Oppure:

Techne Incorporated, University Park Plaza, 743 Alexander Road, Princeton, New Jersey, 08540-6328, USA.

Telefono: (609)452-9275

Toll free: 800-225-9243

Telefax: (609)987-8177

e-mail: salesus@techne.com

Web site: www.techneusa.com

INTRODUCTION

Read the whole of this book before commencing work with the unit.

Blotting techniques in which DNA, RNA or protein are immobilised onto nylon or nitrocellulose filters form a cornerstone of molecular biology. Detection of the desired sequence is usually performed using radioactive probes. Given the expense and hazardous nature of the probes, probing should:

- a) be performed under safe conditions
- b) use minimal volumes (to improve signal and reduce expense)
- c) allow reuse of materials
- d) be simple to use

The Techne Hybrigene meets all these requirements.

Up to five 25 x 25cm filters can be probed in one of the large hybridisation tubes. Hybridisations can be performed in a minimal volume of 5ml, an 80% reduction over conventional systems, reducing expense and improving signal. The operator is protected from irradiation throughout and the apparatus is extremely simple to use. However, the chemical manufacturer's safety precautions must be used when using the Hybrigene.

The Techne Hybrigene also incorporates the following features:

- a) a temperature range of ambient +10°C to 80°C.
- b) it holds 1, 2, 3 or 4 x 80mm diameter hybridisation tubes.
- c) it will hold 3, 6, 9 or 12 x 44mm diameter tubes if you use a three-tube carrier.
- d) it will hold 4, 8, 12 or 16 mini tubes if you use a four-tube carrier.
- e) the Hybrigene comes with adjustable feet to enable accurate levelling.

Initial Checks

When you receive your new Hybrigene remove all the packing. Check the unit for any transit damage. If there is any transit damage at all inform Techne Sales Department or your Techne dealer immediately.

When unpacking the unit, check that the following have been removed from the packing: **a power cable; a tube of silicon grease; a decontamination certificate and a Guarantee Card.**

Check that the unit is the correct rating for your power supply by looking on the Rating Plate which is on the rear of the unit.

Within the guarantee period, shown on the Guarantee Card, we undertake to supply replacements free of charge for parts which may, on examination, prove to be defective, provided that the defect is not the result of misuse, accident or negligence. Any instrument requiring service under this guarantee should be sent to the supplier through whom it was purchased, or, in the case of difficulty, it should be carefully packed in its original packing and consigned, carriage paid, to us. Techne takes no responsibility for returned goods damaged in transit.

Returned goods will not be processed without a Returns Authorisation Number. Call + (44) (0)1223 832401 for a number. On all correspondence, please quote the Serial Number in full and/or the Sales Order Number. **Please write the Returns Number on the outside of any packing.**

Specification


Maximum Number of Hybridisation Tubes	4 (large and small) 16 mini tubes
Temperature Range #	ambient +10° to 80°C (NOTE: minimum temperature must be greater than 10°C above ambient)
Rotation of Hybridisation Tubes	5 rpm to 20rpm and off
4 steps, 5 positions and off	Approx steps of 4 rpm
Adjustable rocking platform	5-20 or 15-60 opm
Temperature stability in chamber	better than 1.0°C
Stability in the tubes	better than +0.1°C
Temperature uniformity in the chamber *	better than +1.5°C
Temperature uniformity in the tubes *	better than +1.0°C
Temperature set point resolution	0.1°C
Absolute accuracy	better than +0.3°C
Heater Power 230V and 115V	750 W
100V	525 W
Temperature selection	Digital
Temperature Display	LED
Temperature Sensor	PRT
External Dimensions	
Width	355mm
Depth	383mm
Height	432mm
Weight	
Net	15 Kg

The Hybridiser includes a preset re-settable overtemperature cutout.

* Temperature measured in the centre of each tube.


Large tubes are fitted with 70°C seals, seals part number 6101488 are required to work to 100°C.

For the Hybridisation Tubes see the list under Accessories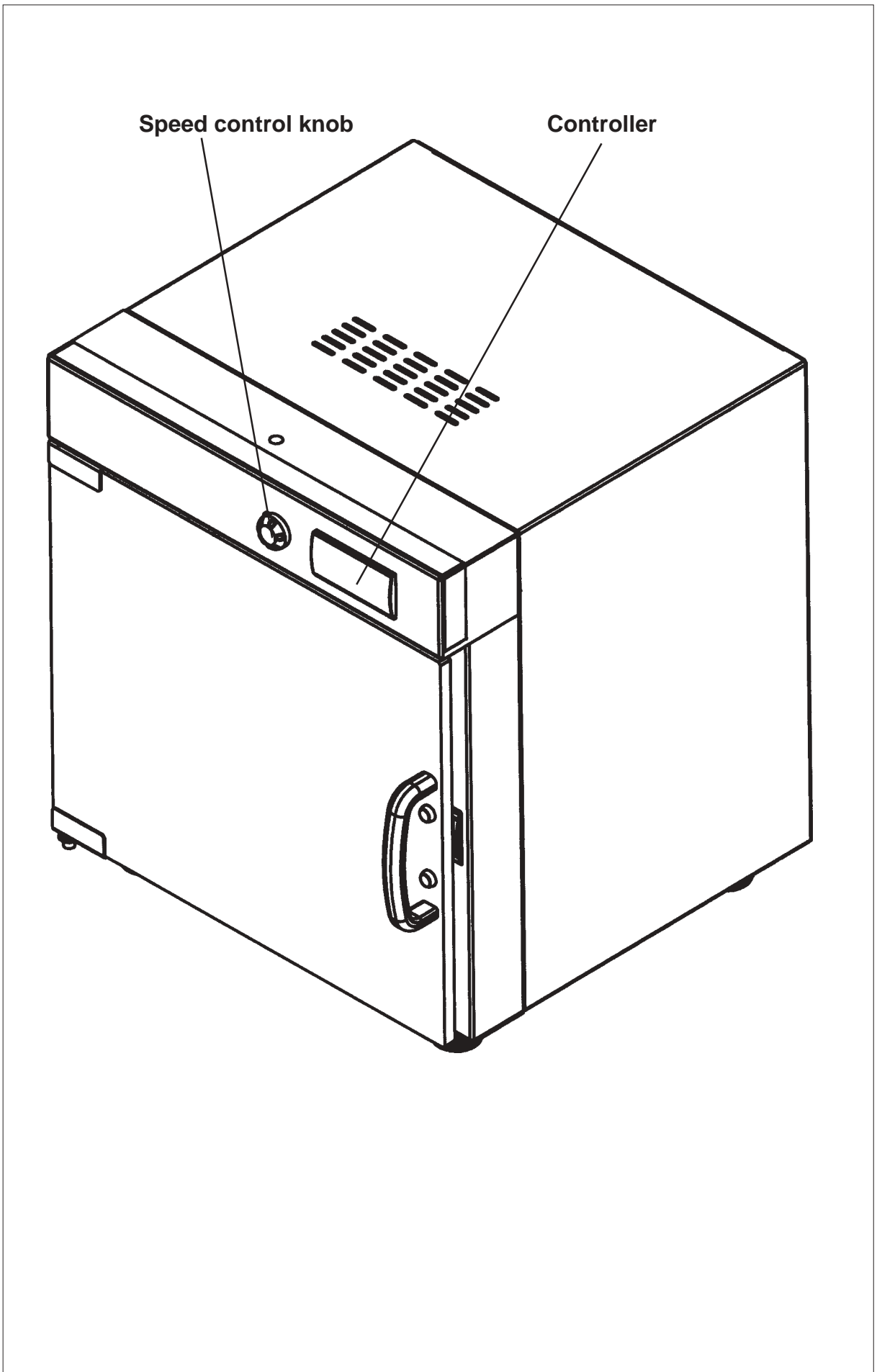

Installation

- 1 Place the unit on a suitable bench, a flat workspace, or in a safety cupboard if required, ensuring that the air inlet and outlet louvres are not obstructed.
- 2 Insert some liquid into a tube and place the tube in the Hybrigene. Adjust the feet by unscrewing them anticlockwise slightly until the unit is level. Check the level by putting the tube in different positions.
- 3 Plug the mains cable into the socket on the rear of the unit.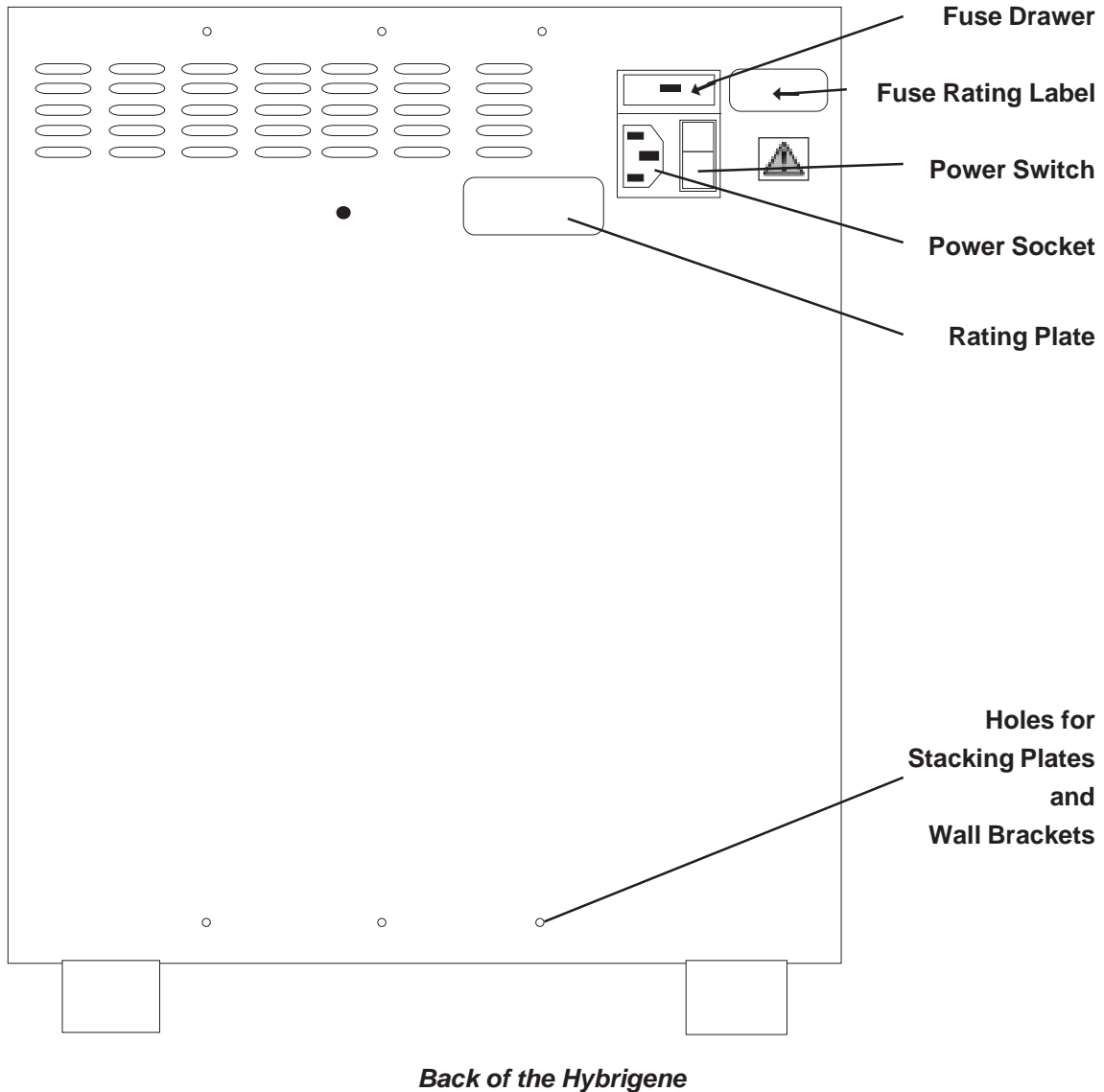

Fitting stacking plates and wall plates on a Hybrigene

A trained technician should carry out this work only after reading the full instructions.

In a **full kit**, check that you have:


- A wall bracket;
- A support bracket;
- 2 stacking plates;
- 15 M4 x 8mm long screws;
- 15 M4 crinkle washers;
- 3 No 8 x 3/4" self tapping screws;
- A thumb screw.

In a **wall bracket kit**, check that you have:


- A wall bracket;
- A support bracket;
- Three M4 x 8mm long screws;
- Three M4 crinkle washers;
- Three No 8 x 3/4" self tapping screws;
- A thumb screw.

In a **stacking kit**, check that you have:

- A stacking plate;
- Six M4 x 8mm long screws;
- Six M4 crinkle washers;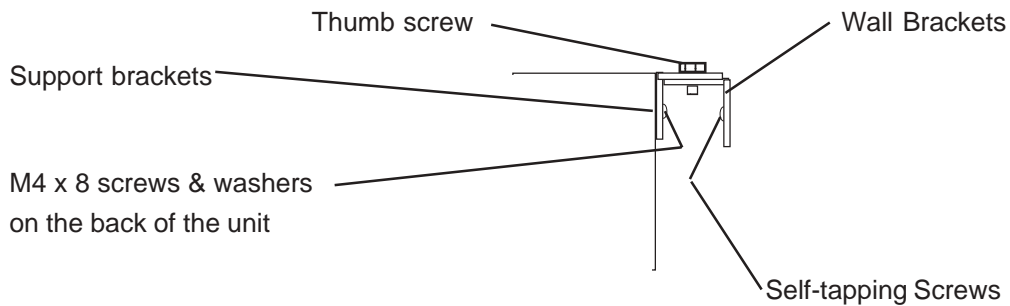

Side of the Top Hybrigene

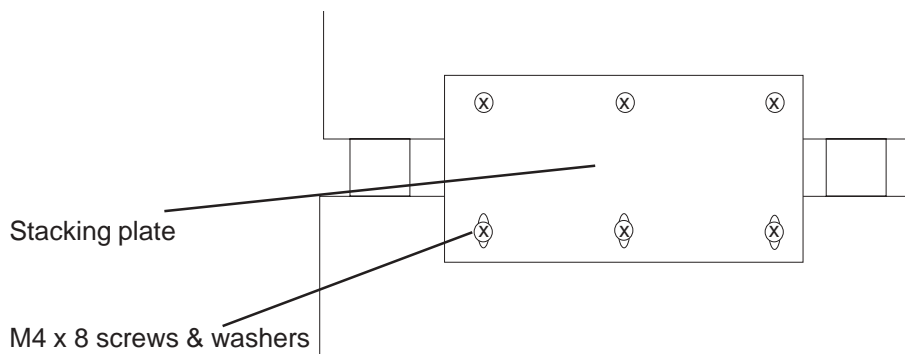Backs of the Hybrigenes

Fixing Procedure

Two Hybrigenes

- Switch the units off and remove the power to them.
- Put two units one on the other.
- The Stacking Plate has three holes and three slots. It has three slots to allow for the adjustment of the feet to get the hybridisers level.
- Use three of the screws to attach the plate to the bottom of the top hybridiser.
- Use the other three screws to attach the plate to the bottom hybridiser.
- Ensure that the screws are tight.

Three Hybrigenes

- If you have three units repeat the process for the top two units. We recommend that you fix the top Hybrigene to the wall for safety.

Wall Fixing

- Use the three screws to attach a support bracket to the top of the top unit.
- Ensure that the screws are tight.
- Put the hybridisers in the position you want to use them.
- Position the wall bracket and mark the position of the holes.
- You should use the appropriate wall fixings for the type of wall you are securing the units against.
- Use the self-tapping screws or the appropriate screws to attach the wall bracket.
- Whatever fixture you use, ensure that the screws are tight.
- Re-position the hybridisers.
- Use the thumb-screw to attach the support bracket to the wall bracket.

Switching On

Once your Hybrigene is connected to a power supply, switch it on using the power switch next to the power inlet socket on the rear panel.


Speed control knob

Controller

Front panel of the Hybrigene

Temperature Control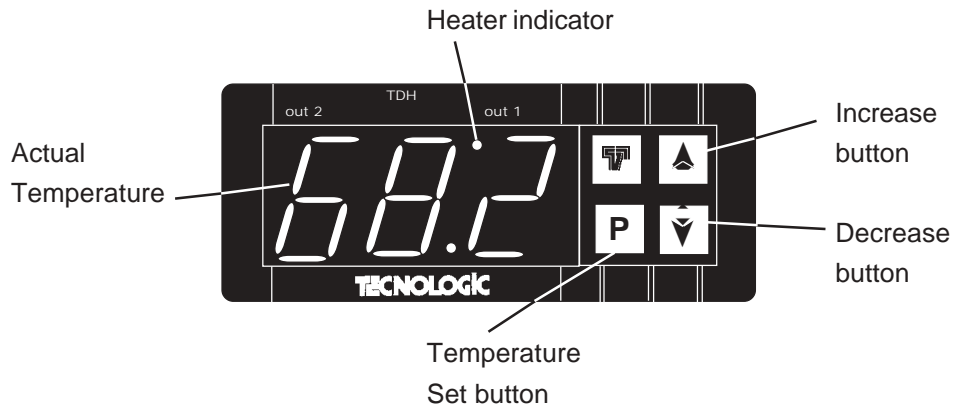

Controller Display

Press the 'P' button on the controller and the set temperature will be displayed.

To increase the set temperature press the up arrow button.

To decrease the set temperature press the down arrow button.

If no key is pressed for 5 seconds the display will return to the measured temperature and the unit will go to the new set temperature.

The Heater Indicator will be lit while the heater is on.

The set temperature can be checked at any time by pressing the 'P' button.

Speed control

Fully anticlockwise the speed control knob switches the rotation of the tubes off

It then increases the speed clockwise and decreases the speed anticlockwise in steps of approximately 4 rpm. Switch the knob until you have the speed you require.

Warning

THE UNIT WORKS UP TO 80°C. This temperature can cause burns. Techne take great efforts to protect operators from hazards, but the following points are brought to your attention:

- 1 Use care and wear protective gloves to remove hybridisation tubes from the cabinet if the unit is at high temperature.

If the protocol allows, it is better to switch off the unit and open the door to allow the unit to cool to below 40°C before attempting to remove the tubes.
- 2 The coloured end cap of the 80mm large tube and the 44mm small tube has two holes so you have to be careful not to spill the probe through these.

One is the hole which allows the air to expand when the tube is heated.

The other is the filling hole which is revealed when the spindle cap is removed.

DO NOT TIP THE TUBE SO THAT THE COLOURED END IS DOWN AS THIS WILL CAUSE A SPILLAGE.
- 3 The unit has been designed to conform to international safety specifications and is fitted with a resettable over-temperature cutout. If, for any reason, the temperature rises above a safe level for the instrument, the overtemperature protection cuts out the heater, but the tubes continue to rotate. Check the reason for the cut out and, if there is no obvious cause, reset the cutout once the unit has cooled. The reset, on the back of the unit, is operated by pressing the centre of the button.
- 4 If the Controller displays 'Err', the thermistor has broken or become disconnected and the unit will cool down automatically.

Servicing by Techne and their Agents

Please note that neither Techne nor any of their agents can accept the return of ANY GLASSWARE from these units.

If the unit is returned it must be thoroughly decontaminated and a CERTIFICATE OF CLEANLINESS returned with the unit.

A copy of the certificate is available from our web-site, www.techne.com, or can be requested from the Service Department if the unit is being returned.

OPERATION

The Tubes


The Hybridisation Tubes must be checked regularly and replaced if any damage occurs. Any damage to the ends of the tubes can cause them to break when the end caps are inserted.

Each tube must be fitted with a radiation shielding end cap if you are using a radioactive probe. These end caps are supplied with all tubes which are supplied with the HYBRIGENE and any other tubes supplied after December 1999. Normal handling procedures must be observed when handling any tube.

Load the membrane and reagents according to which size hybridisation tube you are using.


Large tube (80mm diameter) assembly

Remove the tube end cap which has the holes in it. Place the filters in the tube and replace the end cap. A small quantity of Techne O-ring grease, see the list of accessories, **must** be smeared on the seal to facilitate the insertion of the end cap. Remove the spindle cap and the liquids appropriate to the protocol may then be added through the hole in the spindle using a syringe. Replace the spindle cap and press it onto the spindle so that it clicks into place. Note: the tube will not rotate properly if the cap is not fitted securely.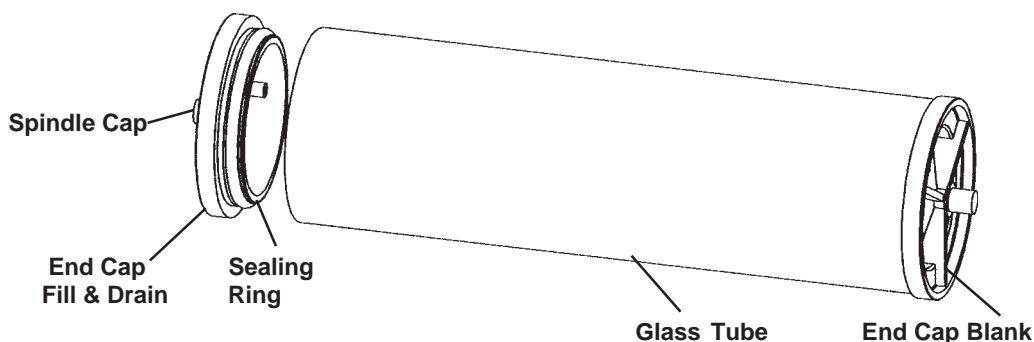

Mini tube (4 x 32mm dia) assembly


Unscrew the end cap of the tube. Place the filters and the liquids appropriate to the protocol in the tube and replace the end cap. Fix the tube into a multi-tube holder, using the springs, and ensure that there are four tubes in the holder. The end caps must be as shown, on alternate ends. The multi-tube holder will not rotate properly if there are not four mini tubes in it.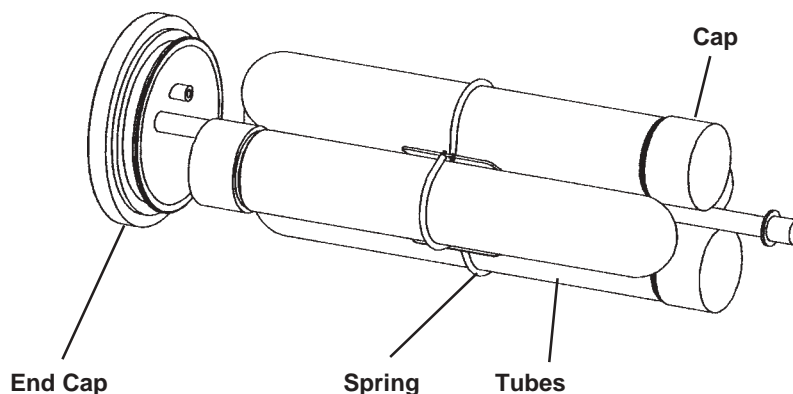

Small Tube (3 x 44mm dia) assembly

Unscrew the tube end cap. Place the filters in the tube and replace the end cap. Remove the spindle cap and the liquids appropriate to the protocol may then be added through the hole in the spindle using a syringe. Replace the spindle cap and press it onto the spindle so that it clicks into place.


The assembly is similar to the four tube assembly but the end caps do not have to be at different ends of the carrier. Fix the tube into a multi-tube holder, using the springs, and ensure that there are three tubes in the holder. The multi-tube holder will not rotate properly if there are not three small tubes in it.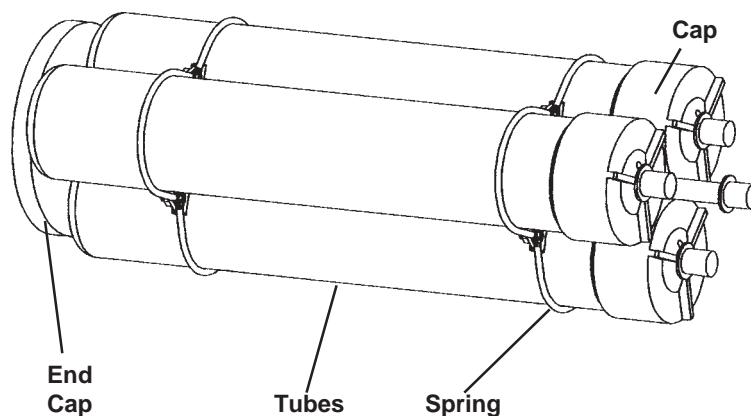

16 x 15ml Tube assembly

Unscrew the end cap from the tube. Place the filters and the liquids appropriate to the protocol in the tube and replace the end cap.

Assemble the tubes into a multi-tube holder, using the spring clips. The multi-tube holder will not rotate properly if it is not balanced properly. Preferably the multi-tube holder should be filled completely even if all the tubes are not being used.

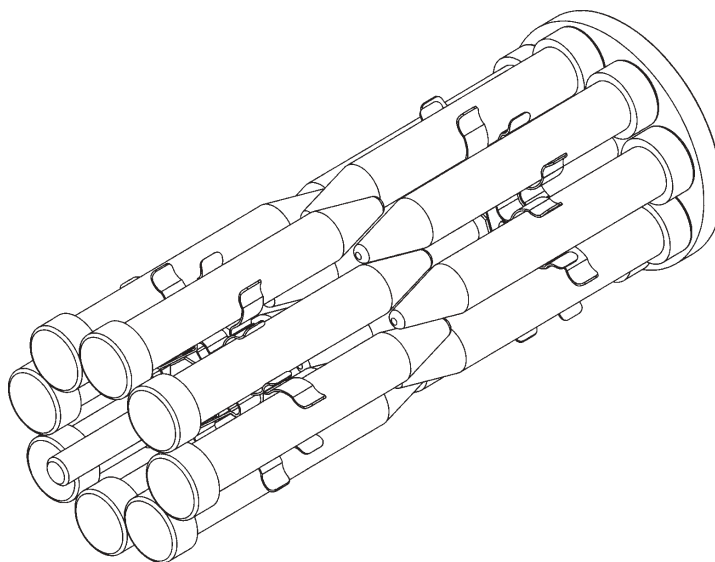The Hybrigene

- 1 Insert the tubes or tube holder into the Hybrigene ensuring that they rotate freely and engage on the drive rollers.
- 2 Close the door.
- 3 Connect the mains electricity supply to the mains socket at the rear of the unit. Switch the power ON. The Controller display on the front panel will light
- 4 Press and hold the top left hand button (the bottom one is inoperative) on the controller and, at the same time, press one of the right hand buttons. The up arrow button will increase the set temperature and the down arrow button will decrease the set temperature.
- 5 The Heater Indicator will be lit while the heater is on.

BASIC PROTOCOL FOR SOUTHERN HYBRIDISATIONS *

INSERTION OF FILTERS

Place a wet or moist filter, DNA side facing inwards, into a hybridisation tube and allow it to adhere to the glass. Large trapped air bubbles should be smoothed out with a glass pipette; smaller air bubbles are acceptable and do not interfere with hybridisation so long as the surface of the filter is covered with hybridisation fluid.

Several filters may be placed in one tube; there is a slight loss of signal, but no increase in background in the lower filters. To improve the signal we recommend the use of separation membranes, part number FMEM2020.

PRE-HYBRIDISATION

Add 5 or 10ml of fluid, according to the number of filters and the size of tube, so that the filters are covered. Allow to pre-hybridise for usual period.

HYBRIDISATION

Drain off the pre-hybridisation fluid by inverting the tube. Add hybridisation liquid with radioactive probe added by using a syringe and needle through the hole in the blue end-cap. The minimum volume for a single filter is 5ml (large tubes).

WASHES

Washing depends on the stringency required.

One method is to wash genomic blots in the following fashion:

- 1) 2 x SSC, 50ml for 10 minutes
- 2) 2 x SSC, 50ml for 10 minutes
- 3) 0.1 x SSC, 0.1% SDS x 50ml for 10 minutes

This gives very clean backgrounds.

NOTE

- 1) All radioactive waste can be collected for disposal; none need be disposed of down the sink.
- 2) Radioactive probes can be collected and reused several times.
- 3) As the protocols for this unit usually involve the use of radioactive materials the appropriate manufacturer's safety precautions must be taken into consideration.

* Courtesy of: Dr M J S Dyer.

When the Protocol is Finished

When the protocol has finished switch off the unit. Open the door and remove the tubes, again being careful to keep the coloured end-cap uppermost. Remember the tubes may be hot. Either wear gloves or, if the protocol permits, allow the tubes to cool before removal.

TECHNICAL INFORMATION

Brief fault finding notes are given in this section.

NOTE THAT THIS EQUIPMENT SHOULD ONLY BE DISMANTLED BY PROPERLY TRAINED PERSONNEL. **REMOVING THE CASE EXPOSES POTENTIALLY LETHAL MAINS VOLTAGES.** THERE ARE NO OPERATOR MAINTAINABLE PARTS WITHIN THIS EQUIPMENT.

Operator Maintenance

General

Should you experience any problems with your Hybrigene which cannot be easily remedied, you should contact your supplier and return the unit if necessary.

Fill in the **Decontamination Certificate** provided with the unit and return it to Techne Service department giving as much information as possible to help diagnose the fault. Please remember to return the unit in its original packing, Techne accept no responsibility for damage to units which are not properly packed: if in doubt, contact your supplier.

Overtemperature Cutout

In the event of no heater power, check the mains plug and lead, then reset the cutout control by switching the unit off and then on again. The cutout should reset itself, if it does not then the unit needs repair.

Repeated operation of the cutout indicates a serious fault: you may need to return your unit to your supplier for repair.

Fuses

If the controller fails to light when power is supplied to the unit the fuses may have blown. Check that there is no external cause such as a faulty plug; check both fuses and replace any faulty fuse with one of the correct value. Fuse values are given on the label next to the mains input socket.

O ring seals on the Hybridisation Tubes

The O ring seals on the 80mm dia Hybridisation Tubes must be checked regularly and replaced before a leak occurs. Each tube is provided with 6 spare O rings and further O rings can be obtained from Techne or your dealer.

Hybridisation Tubes

The Hybridisation Tubes must be checked regularly and replaced if any damage occurs. Any damage to the ends of the tubes can cause them to break when the end caps are inserted. Further Tubes can be obtained from Techne or your dealer.

Cleaning your Hybrigene

Before cleaning your unit ALWAYS disconnect from the power supply and allow to cool.

You can clean the case of the Hybrigene with a cloth dampened in soapy water. No part of the unit should be immersed in solvents. Do not use acetone or abrasive cleaners.

DO NOT USE ETHANOL OR METHYLATED SPIRITS ON THE ACRYLIC DOOR; these may make the acrylic go cloudy.

If You Spill the Fluid

Because of the nature of operation of the Hybrigene, the complete inside must be cleaned if a spillage does occur. This is to remove any airborne particles which may have been blown around the system and to guarantee full operator safety.

Switch the Hybrigene off and disconnect it from the power supply. Follow this sequence of operations when the Hybrigene has been allowed to cool below at least 50°C :

- a) Remove all glassware from the Hybrigene.
- b) Remove the Plastic drip tray.
- c) Clean all surfaces of the inner chamber using a suitable cleaning and decontaminating solution.

Be careful of the drive shafts.

Cleaning and decontamination materials which Techne recommend include:

Decon 90

Neutracon

or mild detergent.

Accessories and Spares

The following may be purchased from Techne or their dealers. All Techne tubes are made from Borosilicate glass. These tubes are the preferred tubes (they also fit the HB-3D):

FHB16	80mm external diameter hybridisation tube assembly complete with sealing rings and end caps
FHB15	80mm external diameter tube only
6101082	End cap with holes (blue) less sealing ring
6101387	Spindle cap (for use with 6101082).
6101083	End cap without holes (natural) less sealing ring
6101090	Sealing ring (for 70°C operation)
6101351	Techne O-ring Grease (tube)
FHB41	4 x 31mm external diameter tubes only
FHB61	4 tube holder
6102375	End cap for the mini tube
6102376	Seal for the mini tube
6101387	Spindle cap
6101828	Spring for the multi-tube holder
FHB62	8 tube carrier (Falcon disposable tubes)
FHB72	3 tube carrier assembly complete with seals and end caps
FHB71	Tube 44mm diameter
FHB70	3 tube carrier
6101387	Spindle cap
6007625	Seal (for use with 44mm external diameter tubes).
6102193	Screw on end cap, less seal

FHTRACK	Hybridisation Tube Rack
FMEM2020	Separation Membrane 20cm x 20cm
FHB4ST1	Rocking Platform 20/40rpm
FHB4ST2	Rocking Platform 20/60rpm
FHB4SH1	Shelf

Contact Techne Sales or your dealer for further details.