

Media Release

DataWay PC Software

An exciting new software package to accompany Jenway instruments

- ◆ Automatic detection and set-up
- ◆ Log data in tabular or graphical form
- ◆ Direct export to Excel spreadsheets
- ◆ Logs all available data from instrument mode with date & time
- ◆ Data table, graphical or split screen display
- ◆ Store, print or transfer results
- ◆ Easy to use
- ◆ On-Screen prompts and detailed Help files
- ◆ Low cost

Jenway have launched DataWay, a new PC software program for data acquisition for Jenway spectrophotometers and electrochemistry products. DataWay is an expandable, PC based, data acquisition software that is compatible with Windows 98, 2000, ME, and XP.

DataWay automatically recognizes the instrument connected to the PC. No manual settings are required. A "quick start" option enables specific instruments to be selected, speeding up the search and detect process.

Data can be downloaded from the instrument to the PC manually or on a time (data logging) basis. Automatically or manually logged data can be displayed in both tabular and graphical form and a single click exports the data directly to an Excel® spreadsheet. The data is shown along with date and time with an option to display either the data or the graph, or split screen so that both can be viewed together.

Data stored on the PC can be neatly organized in relevant folders accessed through the familiar explorer screens. Expanded print options allow the printing of graph and table on separate sheets or together with either the graph or table on top.

DataWay is intuitive and easy to use, with on-screen prompts and detailed Help files. It can acquire data from all modes of the instrument. Familiar Toolbar icons give quick access to most functions and the "tabbed page" format clearly identifies main working areas. DataWay meets all of your lab needs: a single click transporting data to a spreadsheet for further manipulation, timed data logging and expanded print options is now possible.

DataWay PC software supports the following Jenway instruments: 3510, 3520, 4520, 3540, 6400, 6405, 6500, 6505, and Genova

ORDER CODE: 050-501

DataWay software on CD-ROM includes spectrophotometer interface cable (013-210) and electrochemistry interface cable (013-203)

CONTACT:

Techne Incorporated

3 Terri Lane, Suite 10 Burlington, NJ 08016,
Phone: 800-225-9243
www.techneusa.com

