TECHNE

TC-412

OPERATOR'S MANUAL

Barloworld
Scientific

Issue 1
2/07

CONTENTS

	PAGE
INTRODUCTION	1
SAFETY AND INSTALLATION	2
SÉCURITÉ ET CONSIGNES D'INSTALLATION	5
SICHERHEITS- UND INSTALLATIONSINFORMATIONEN	8
INFORMACIÓN DE SEGURIDAD E INSTALACIÓN	11
INFORMAZIONI SULLA SICUREZZA E L'INSTALLAZIONE	14
BEFORE SWITCHING ON	17
CONTACT INFORMATION	18
TC-412 SPECIFICATION	19
USER OPERATION AND CONSUMABLES	21
Heated lid	21
Tubes or reaction vessels	22
Switching on	22
FRONT PANEL CONTROLS	23
The LCD display	23
Indicators	23
The keys	24
PROGRAMMING THE TC-412	25
Programming Options	26
Option 1: Creating programs using templates	26
Option 2: Creating a new program	30
Example programs	31
Programming	31
To program increment or decrement temperature or time.	34
To run a program	38
INFORMATION	40
CONFIGURATION	40
RUNNING THE TC-3000 FROM A PC	43
Connecting one unit to the computer using an RS232 cable	43
Connecting several units to the computer using the daisy chain cables	44
Upgrading the software in your TC-412	45
USER MAINTENANCE	45
Cleaning your TC-412	45
Block removal	46
FREQUENTLY ASKED QUESTIONS	46
ADDITIONAL INFORMATION	47
Fault Finding	47
Fuses	47
The heated lid over-temperature cut-out	48
Insulation testing	48
INTERCHANGEABLE BLOCKS	48
ACCESSORIES	49
REPLACEMENT PARTS	49

INTRODUCTION

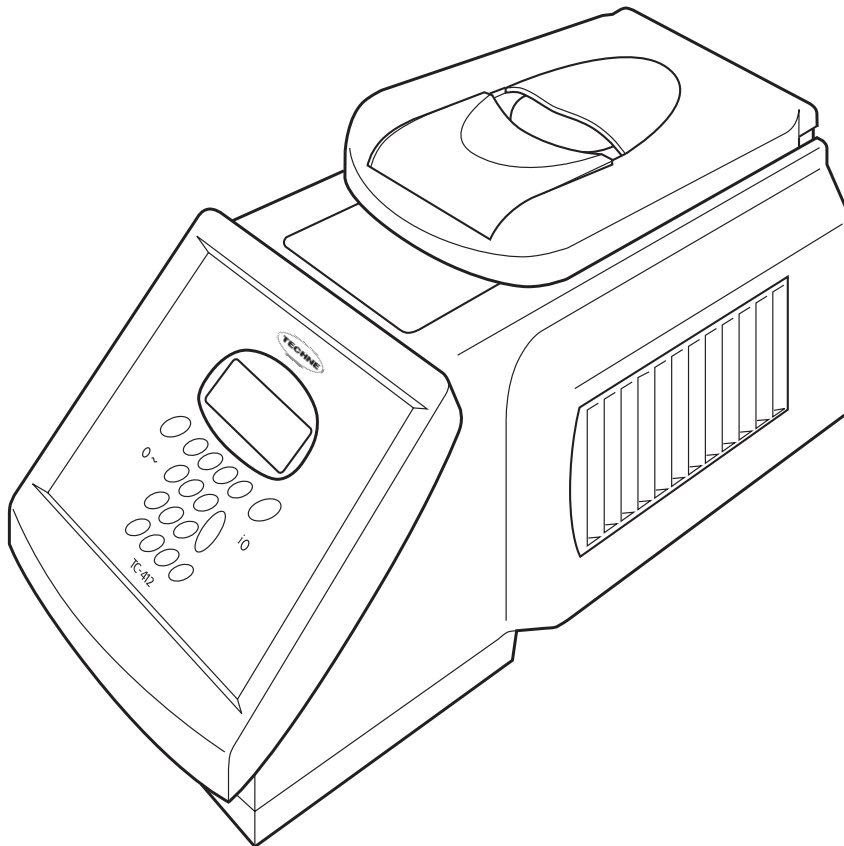
The TC-412 provides the researcher with the means of accurately controlling the temperature profile of samples. It has many scientific applications, including DNA amplification¹ and sequencing. The TC-412 can cycle samples between 4°C and 99°C.

The TC-412 is programmed by means of an integral keypad and LCD display. A program, which can be recalled from memory, consists of:

- a series of specified temperatures in °C
- the times for which each specified temperature will be held (Hold Times)
- the desired heating or cooling rates, in °C/sec, between each specified temperature
- optional instructions as to whether the times and/or the temperatures are to increase or decrease when a cycle is repeated.

The memory can store up to 80 programs.

Before using the TC-412 make sure you have read this manual carefully. If there is any doubt relating to the proper use of this equipment, the staff at Barloworld Scientific Ltd or your supplier will be happy to assist you.

¹Aspects of the PCR process are claimed in USA Patent Nos. 4,683,195; 4,683,202, and 4,965,188. Use of the TC-412 in such processes does not convey a licence to practice the processes themselves.

SAFETY AND INSTALLATION

OPERATOR SAFETY

Please read this manual carefully before using the thermal cycler. If the equipment is not used in the manner described in this manual the protection provided by the equipment might be impaired.

All Barloworld Scientific Ltd instruments are designed to conform to international safety requirements and are fitted with an over-temperature cut-out. If a safety problem should be encountered then switch off the unit at the mains socket and remove the plug from the electricity supply.

WARNING

HIGH TEMPERATURES ARE DANGEROUS as they can cause serious burns to operators and ignite combustible material. Users should be aware of the following potential hazards:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS
- DO NOT use combustible substances near hot objects
- DO NOT operate the instrument in the vicinity of inflammable liquids or gases
- DO NOT place any liquid directly into the instrument, always use the appropriate plastic consumables

After heating samples, remember that parts of the instrument, namely the lid and block and any associated accessories may be very hot.

UNPACKING

When unpacking please ensure that the following have been removed from the packaging:

- Thermal cycler
- Mains cable
- Guarantee card
- Decontamination certificate

The user is advised to keep the original packaging in case the instrument ever needs to be returned for service or repair. Barloworld Scientific Ltd accepts no responsibility for damage incurred unless the unit is correctly packed and transported in its original packaging.

Please note that the screen of the thermal cycler is easily damaged by sharp objects such as pens, pencils and fingernails. This type of damage will be considered as misuse and invalidate the guarantee for this component.

INSTALLATION

The instrument should be carried using both hands. Never move or carry the instrument when in use or connected to the mains electricity supply.

1. All Barloworld Scientific Ltd instruments are supplied with a power cable; this may be integral or plug-in.
2. Before connecting the instrument to the mains electricity supply, check the voltage against the

rating plate (located on the back of the unit). Ensure that the voltage selector switch (located above the On/Off switch) is set to the appropriate voltage for the local supply. Please note that the unit must be earthed to ensure proper electrical safety. Connect the mains cable to a suitable plug according to the table below.

Connections	220V-240V, 50/60Hz Supply	110V-120V, 50/60Hz Supply
Live	brown	black
Neutral	blue	white
Earth	green/yellow	green

3. UK ONLY: The fused plug supplied with the mains cable is fitted with a 10 Amp fuse to protect the instrument and the user.
4. Units showing 230V, 50/60Hz on the rating plate also operate between 198 and 252V; units set to 115V, 50/60Hz also operate between 110 and 130V and units marked with 100V will operate between 90 and 110V. In all cases the heating rate will degrade by approximately 8% at the extremes of the voltage range.
5. Plug the mains cable into the socket on the back of the instrument.
6. Place the unit on a suitable flat bench or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction. Position the instrument with a minimum distance all round of 200mm from walls or other items and between each unit if multiple units are being used to permit efficient air flow for each instrument.
7. Switch on the instrument:
 - mains switch On
 - mains switch Off

Replacement cable

Should the mains lead need replacement, a cable of 1mm² of harmonized code H05VV-F connected to an IEC320 plug should be used. **IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN.**

WORKING CONDITIONS

The thermal cyclers are designed to operate under the following conditions:

- Indoor use
- Ambient temperature range +5°C to +40°C
- Altitude to 2000m
- Mains supply fluctuations not exceeding 10%
- Over voltage category II IEC 60364-4-443
- Pollution degree 2

Note: The control specifications are quoted at an ambient temperature of 20°C. The specification may deteriorate outside an ambient temperature of 10°C to 30°C.

The instrument has been tested for radio frequency interference and is certified under EN61326.

GUARANTEE

The instrument is guaranteed against any defect in material or workmanship for the period specified on the enclosed guarantee card. This period is effective from the date of purchase; within this period all defective parts will be replaced free of charge provided that the defect is not the result of an accident, misuse or negligence. The screen of the thermal cycler can be damaged by sharp objects such as pens, pencils and fingernails. Damage of this sort will be considered misuse and will invalidate the guarantee for this component.

Servicing under this guarantee should be obtained from the supplier of the instrument.

This manual has been prepared for the convenience of Techne's customers and nothing in this manual shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the unit or components.

Notwithstanding the description and specification(s) of the instruments contained in the operator's manual, Techne reserves the right to make such changes as it sees fit to the instruments or to any of the components.

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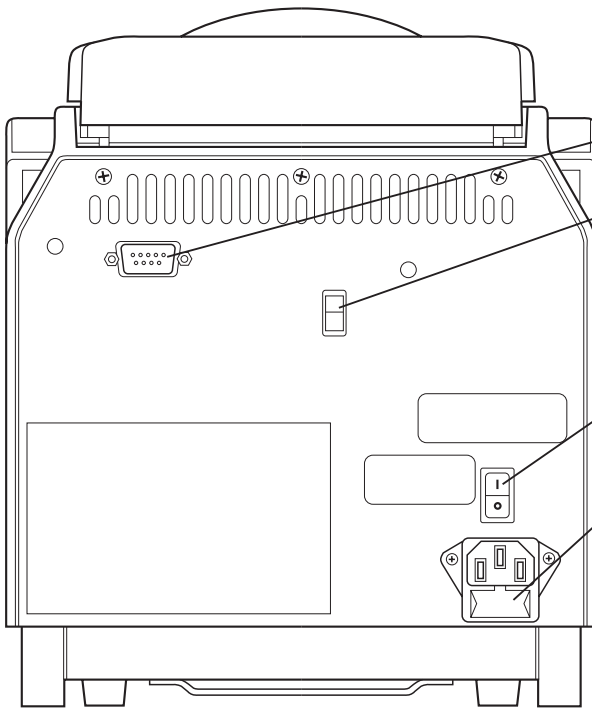
BEFORE SWITCHING ON

Serial port

Ensure that the voltage selector switch is set to the appropriate voltage for your supply;

On/Off switch

Ensure that the appropriate fuses are fitted for your supply:

- For 230V set the voltage selector switch to 230V and fit T6.3A fuses.
- For 115V set the switch to 115V and fit T8A fuses.
- For 100V the switch will be set to 115V; T10A fuses will be fitted.

CONTACT INFORMATION

For technical, sales or servicing information, contact your local Barloworld Scientific Ltd dealer or:

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TC-412 SPECIFICATION

Temperature

Temperature range	4°C to 99°C
Temperature set point precision	0.1°C
Block uniformity (over full range)	± 0.6°C
Temperature accuracy	± 1.0°C
End of program cooling below ambient	Yes

Heating/Cooling Rate

Heating rate, between 55°C and 90°C on a 50°C to 95°C step:	
96 x 0.2 ml micro tubes	2.6°C/sec
60 x 0.5 ml micro tubes	2.5°C/sec
Cooling rate, between 90°C and 55°C on a 95°C to 50°C step:	
96 x 0.2 ml micro tubes	1.3°C/sec
60 x 0.5 ml micro tubes	1.6°C/sec
The programmable ramp rate is settable to a maximum of 2.0°C/sec. Any value above 2.0°C/sec will default back to 2.0 unless set for max. If the unit is set to max, the heat up rate will be 2.6°C/sec.	

Heated Lid

Heated lid enable/disable	Yes
Selectable heated lid temperature	100°C to 115°C
Over-temperature cut-out	Fixed at 145°C
Temperature sensor	Thermistor
Heater type	Wire wound
Heater power	132W
Warm up time	3 min approximately (ambient to 105°C)

The heated lid will only come on if the set temperature is above 35°C

Programming

Program naming using alpha keys	Yes
Program password protection	Yes
Number of programs (3 step)	80
Maximum number of steps per stage	35
Maximum number of stages per program	25
Minimum hold time	1 sec
Maximum hold time	18 hrs
Incremental/decremental temperature	Yes
Incremental/decremental step hold time	Yes
Programmable ramp rate (heat or cool)	Variable in 0.1°C/sec steps
Programmable ramp rate resolution	0.1°C

Running Programs

Pause button	Yes
Stop button	Yes
End of program alarm	Yes
Auto restart on power fail	Yes

Interconnectivity

Connection to PC control program	Yes
Software updates from Techne web-site	Yes

Dimensions

Height	260mm
Width	220mm
Length	420mm

Serial Port

This socket is provided for RS232 PC link to run "Gensoft" the PC control software.

Power

Power consumption	620W
-------------------	------

The units are voltage selectable by a slide switch. When the unit is set to 115V it will work at any voltage between 110V and 130V. When the unit is set to 230V it will work at any voltage between 198V and 252V.

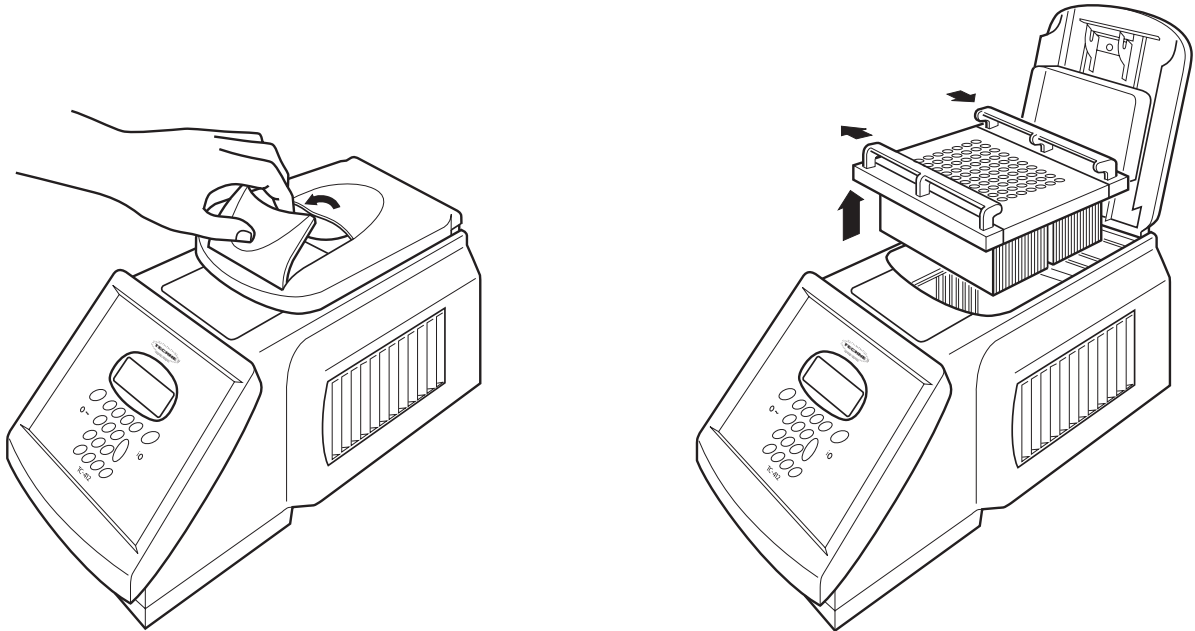
Units marked 100V will work at any voltage between 90 and 110V.

The performance will vary and will not necessarily meet the above typical specification at the extremes of voltage.

USER OPERATION AND CONSUMABLES


HEATED LID

To release the heated lid, lift the back of the handle as shown and use it to lift the lid. To close the lid, lower the lid onto the block containing the samples and press the handle down to shut.

The silicone rubber mat supplied with the unit (Part No. 6104978) is for use with microplates when sealing using a plastic adhesive film. The mat should be placed on top of the sealing film to ensure good thermal contact with the heated lid.

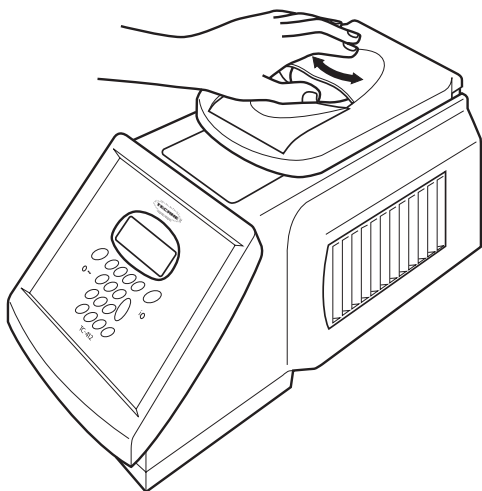The heated lid has a rotating knob to adjust the lid pressure on the samples, allowing for a variety of consumables to be used. To adjust the pressure:

1. Place the samples in the block and place the lid in the down position.
2. Adjust the orange rotating knob anticlockwise to give the least pressure then close the lid fully.
3. Gently rotate the knob clockwise until you can feel pressure beginning to be applied.
4. Rotate the knob a further quarter of a turn and this will give the correct pressure.

Do not over-tighten the knob as this will cause deformation of the silicone rubber membrane (where used) and reduce the effectiveness of the lid. Once the pressure has been set it does not need to be adjusted unless a different block or consumable is used.

TUBES OR REACTION VESSELS

Techne does not recommend any specific tube or reaction vessel other than those described in this manual. We recommend using reaction volumes between 20 and 200 μ l. The tubes must withstand a pressure of 1 atmosphere at 100°C. Any vessel must be able to withstand the temperatures you are using without any danger of them deforming to the point where they fracture.

To test your tubes, put 25 μ l of water into each of 5 tubes and subject them to a typical thermal cycling protocol. At the end of the program, measure the volume remaining using a micropipette. A loss of more than 1 or 2 μ l indicates a vapour leak.

The amount of volume loss you observe and the change in reactant concentrations you can tolerate determine the minimum volume that can be used. Typical volume losses of 1 μ l in 30 cycles allow the use of samples of 20 μ l or less.

During the final cool-down, a ring of condensation may form above the liquid level but below the top of the sample block. This is normal and does not effect the reaction.

SWITCHING ON

```

TC-412
Version          xx.xx
Block           96x0.2ml
                60x0.5ml
                384
                Flat
  
```

When you switch on the unit a screen similar to the one shown here will appear briefly. This indicates the software version and block type installed in the unit.

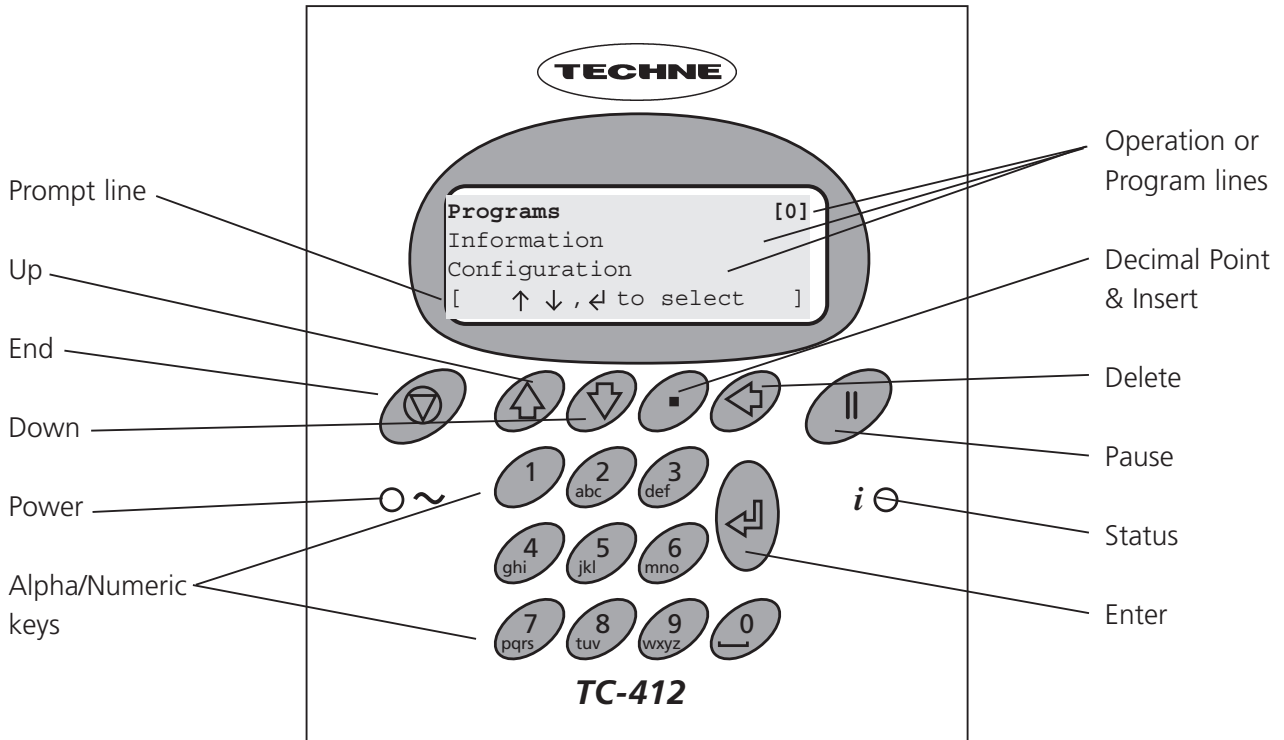
```

Programs [0]
Information
Configuration
[  ↑ ↓ , ← to select  ]
  
```

When the initial screen has disappeared the screen shown here will automatically appear. There are three options:

1. 'Programs' for creating and running programs;
2. 'Information' for obtaining information about the unit e.g. serial number;
3. 'Configuration' for setting up program defaults and settings.

Each of these is explained in detail in the following sections.

FRONT PANEL CONTROLS

THE LCD DISPLAY

```

Programs [0]
Information
Configuration
[ ↑ ↓ , ← to select ]
    
```

The TC-412 has a four line LCD display. The top three lines are information or program lines, while the bottom line is a prompt line. The active line flashes on screen and is shown in bold in this manual.

```

----- Stage 1 -----
Number cycles          10
Step 94.0°C 00m30 max
[ '.' to insert step ]
    
```

The bottom line may change or flash between two or three different prompts depending upon which program line is active. An example of a prompt line which flashes between three different prompts is shown here.

```

[ ↑ ↓ , ← to edit ]
[ ← to delete step ]
    
```

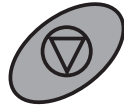
INDICATORS

○ ~ Power: This indicator is lit when there is power to the TC-412.

i ○ Status: This indicator is lit when the TC-412 is running a program. It is off when the unit is stopped. It will flash and beep slowly when the unit is paused. It will flash after a five second quick sequence of beeps when the unit reaches its final hold temperature.

THE KEYS

The End, Finish or Exit key

The "End" or "Finish" key is used to finish editing programs or to exit a program. It is also used to exit a sequence of key operations.

The Up Arrow key

The "Up Arrow" key is used for scrolling up through the lines on the screen. It is also used to increase certain fields when editing.

The Down Arrow key

The "Down Arrow" key is used for scrolling down through the lines on the screen. It is also used to decrease certain fields when editing.

The Decimal Point and Insert key

The "Decimal Point" key is used when editing numeric fields. It is also used as an Insert key during program editing to insert additional steps and other field sets.

The Delete key

The "Delete" key is used during program editing to delete steps and stages.

The Pause key

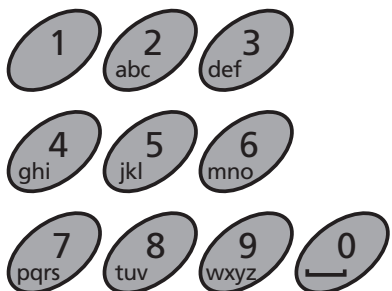
The "Pause" key is used during program editing to insert a pause after a stage or to insert increment and decrement times and temperatures. It is also used to pause a program when it is running.

The Enter or Accept key

The "Enter" key is used to enter an operation or the edit mode. It is also used to 'Accept' a field or program change.

The Alpha/Numeric keys

The "Alpha/Numeric" keys are used to enter numbers and letters during programming. If it is a numeric field then numbers will be inserted; if it is an alpha field then letters will be inserted. Each key will cycle through the letters shown on the key (0 will insert a space). To insert a number, press and hold the appropriate key.

PROGRAMMING THE TC-412

The table below illustrates the symbols which appear on the LCD display and the relevant keypad button to which they refer.

	In the prompt line is equivalent to the "End" key.	
	On a program line means the program is locked; it can be copied and then edited but not edited directly.	
	Up arrow	
	Down arrow	
	Enter	
	Pause	

The following terms are used when programming the TC-412:

- Stage** A 'Stage' is the name given to the main repeated part of a program containing the 'Steps' to be repeated. A program can contain a maximum of 25 stages although typically only one or two are used.
- Cycle** A 'Cycle' is the number of repetitions of a 'Stage'.
- Step** A 'Step' is a user defined program line containing a temperature, a hold time and ramp rate. Typically three steps are created within one 'Stage'.

```

Programs [0]
Information
Configuration
[   ↑ ↓ , ← to select ]
    
```

To enter the **programming** section of the TC-412 software, highlight 'Programs' using the "Up Arrow" or "Down Arrow" keys ↑ ↓

Press "Enter" ←

```

* * New Program * * [ ]
2 STEP TEMPLATE    501  [ ]
3 STEP TEMPLATE    502  [ ]
[   ↑ ↓ , ← to select ]
    
```

This screen lists all the programs saved on the unit. Use the "Up Arrow" and "Down Arrow" keys to scroll through the list. ↑ ↓

The number next to the program name is the unique "Speed dial" number allocated to the program for rapid retrieval (see below).

PROGRAMMING OPTIONS

There are two options available for creating programs on the TC-412:

1. Copy and edit one of the pre-written templates.
2. Create a new program from the defaults.

OPTION 1: CREATING PROGRAMS USING TEMPLATES

There are two pre-written basic PCR templates: a two-step and a three-step template. These can be copied and modified as required.

```

2 STEP TEMPLATE
Heated Lid          105°C
Preheat Lid         on
Pause               off
-----
In denat            94°C 5m00
Hot Start           off
----- Stage 1-----
Number cycles       30
Step 94.0°C 0m30 max
Step 68.0°C 1m00 max
-----
Final Extn 72.0°C 5m00
Final hold          10.0°C
    
```

```

3 STEP TEMPLATE
Heated Lid          105°C
Preheat Lid         on
Pause               off
-----
In denat            94.°C 5m00
Hot Start           off
----- Stage 1-----
Number cycles       30
Step 94.0°C 0m30 max
Step 55.0°C 0m30 max
Step 72.0°C 0m30 max
-----
Final Extn 72.0°C 5m00
Final hold          10.0°C
    
```

Other templates available include Ice bucket (incubation at 4°C), Ligation (incubation at 15°C) and an RT-PCR template.

```

* * * New Program * * *
2 STEP TEMPLATE    501  [F]
3 STEP TEMPLATE    502  [F]
[  ↑ ↓ , ← to select  ]
    
```

Use the "Down Arrow" key ↓ to select an existing 'Template' program from the menu, e.g. '2 Step Template' or '3 Step Template'.

Press the "Enter" key ↵ to accept it.

```

2 STEP TEMPLATE [501]
Run program
View program
[  ↑ ↓ , ← to select  ]
    
```

The template is already defined and can be run as it is or easily copied, modified and saved under a new name.

```

Run program
View program
Copy program
[  ↑ ↓ , ← to select  ]
    
```

Use the "Down Arrow" key ↓ to select 'Copy Program'

Press the "Enter" key ↵ to accept it.

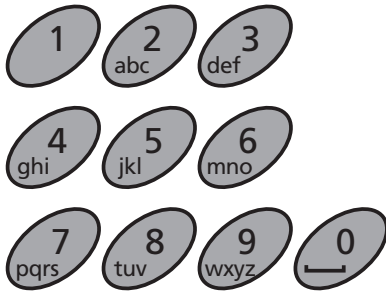
```

UNNAMED #
Heated Lid          105°C
Preheat Lid         on
[   ↑ ↓ , ← to edit   ]
    
```

To give a program a name

```

■ UNNAMED #
Heated Lid          105°C
Preheat Lid         on
[   ↑ ↓ , ← to edit   ]
    
```

```

DONNA
Heated Lid          105°C
Preheat Lid         on
[   ↑ ↓ , ← to edit   ]
[   ▽ to save        ]
    
```

To save the program

```

DONNA
Heated Lid          105°C
Preheat Lid         on
[   SAVE?           ]
[   ← YES or ▽ NO   ]
    
```

```

DONNA [7]
Run Program
View Program
[   ↑ ↓ , ← to select ]
    
```

A new name can now be given to the copied program (see below).

To type in a name for the program, press "Enter" ← and a cursor will appear in the position shown.

Using the keypad, type in the name you wish to give the program.

The letters will scroll round with successive presses; for example the "2" key will give 'A' with one press; 'B' with two presses; 'C' with three presses and '2' with four presses. Five presses will again give 'A'.

When the name is complete, press "Enter" ←

There are now two options:

1. Save the program and run it as it is or
2. Edit the program then save and run it.

A program can only be run once it has been saved. Press the "End" ▽ key.

The prompt line will flash between 'SAVE?' and 'YES or NO'.

Press "Enter" ← for 'Yes' to save the changes you have made.

If you press "End" ▽ for 'No' then all the changes you have made will be lost.

The program is now ready to run and will be allocated with a unique 'Speed dial' number which can be used to rapidly locate the program from the main start-up screen.

The "Up Arrow" or "Down Arrow" keys ↑ ↓ can be used to scroll through the remaining list of options.

To edit a field or step in an existing program

```

Programs [0]
Information
Configuration
[ ↑ ↓ , ← to select ]
    
```

```

DAVID 8 [D]
DONNA 7
[ ↑ ↓ , ← to select ]
    
```

When the unit is switched on the display will show this screen.

If known, type in the 'Speed dial' number for the required program and press "Enter" ← This will take you directly to the required program.

If you do not know the 'Speed dial' number, press "Enter" ← to access the list of stored programs.

Either:

1. Scroll to the required program using the "Up Arrow" or "Down Arrow" keys ↑ ↓ or
2. Type the first letter of the program name in the 'New program' field; this will then list all the program names beginning with that letter. Use the "Up Arrow" or "Down Arrow" keys ↑ ↓ to select the required program.

Press "Enter" to accept ←

```

DONNA [7]
Run program
View program
Edit program
Copy program
Delete program
Set password
[ ↑ ↓ , ← to select ]
    
```

The full set of options for a program is shown here; these are shown on the LCD screen three lines at a time.

Use the "Down Arrow" and/or the "Up Arrow" keys ↑ ↓ to select the option you require.

Press the "Enter" key ← to select the option.

Edit a selected field or step

```

DONNA
Heated Lid 105°C
Preheat Lid on
[ ↑ ↓ , ← to edit ]
[ ▽ to save ]
    
```

Select 'Edit program' and a screen similar to this will appear.

Use the "Down Arrow" and/or the "Up Arrow" keys ↑ ↓ to select the field or step you want to edit.

You can now edit the active field or step by first pressing the "Enter" key ←

```

DONNA
Heated Lid      █ 105°C
Preheat Lid      on
[   edit temperature   ]
    
```

[← to disable]

```

DONNA
Heated Lid      105°C
Preheat Lid      █ on
[   ↑ ↓ to change   ]
    
```

[turn on/off]

```

Pause          off
-----
In denat  █ 94.0°C   00m30
[   edit temperature   ]
    
```

[← to disable]

```

Pause          off
-----
In denat  94.0°C   █ 00m30
[   edit time   ]
    
```

[.for hrs/mins]

```

-----
Fin extn 72.0°C   5m00
Fin hold      10.0°C
[   ↑ ↓, ← to edit   ]
    
```

[▽ to save]

To edit a temperature, type in the required temperature. In the example here, the heated lid temperature is being edited.

Press "Enter" ↵ to accept.

Use the "Down Arrow" and/or the "Up Arrow" keys ↑ ↓ to select the next field or step you wish to edit.

To edit a field with on/off settings, first select the field and press "Enter" ↵

Use the "Up Arrow" and "Down Arrow" keys to toggle ↑ ↓ between the settings.

Press "Enter" ↵ to accept.

To edit a step with both a temperature and a time, first press "Enter" ↵ and type in the new temperature. Press "Enter" ↵ again. Now type in the new time. The default is minutes.

1. For whole minutes, type the number required then "Enter" ↵
2. To toggle between minutes and hours, use the "Decimal Point" key.
3. For values less than 1 minute, first press the "Decimal Point" key followed by the number of seconds required.
4. For minutes and seconds, type the number of minutes followed by the "Decimal Point" key then the number of seconds.

Once the values have been added, press "Enter" ↵ to accept.

Once you have finished editing and you are happy with the changes, press the "End" key ▽ followed by the "Enter" key ↵ to save.

If you press the "End" key ▽ a second time for 'No' then all the changes you have made will be lost.

To copy a selected program

```

                UNNAMED #
Heated Lid           105°C
Preheat Lid           on
[   ↑ ↓ , ← to edit   ]
[           ▽ to save  ]
    
```

If you select 'Copy program' a screen similar to this will appear and the software will automatically give the program the next number until it is named.

Press "Enter" **↵** and type in a new name. You will then be able to edit the program as previously described.

To delete a selected program

```

                DONNA
                delete program
                SURE?
[   ← Yes or ▽ No   ]
    
```

If you select 'Delete program' a screen similar to this will appear.

Press "Enter" **↵** to delete the program or "End" **▽** to return to the program screen.

To set a password for a selected program

```

                DONNA
New passwd   █
[   ← to accept   ]
[           ▽ to exit   ]
    
```

If you want to password protect the program so that other users cannot edit it, select 'Set password'. A screen similar to this will appear. Type in the password you wish to give it and press "Enter" **↵**. The program list will then display the key symbol **█** next to the name to show that it is locked. Other users will still be able to copy/view your program but not edit it.

OPTION 2: CREATING A NEW PROGRAM

A second option is to create a completely new program without using templates.

```

                UNNAMED #
Heated Lid           105°C
Preheat Lid           on
Pause                off
-----
In denat      94.°C      5m00
Hot Start                off
-----
Final Extn    72.0°C      5m00
Final hold                10.0°C
    
```

A number of typical values are pre-entered including:

- Heated lid temperature
- Preheat lid
- Initial denaturation
- Hot start
- Final extension
- Final hold

These values can be edited and the additional stages and steps required for cycling can be added by following the instructions detailed on the following pages.

EXAMPLE PROGRAMS

A couple of example thermal cycling programs are shown below.

Typical 3-Step PCR

```

Heated Lid          105°C
Preheat Lid         on
Pause               off
-----
In denat 94.0°C    5min
Hot Start           off
----- Stage 1 -----
Number cycles       30
Step 94.0°C 0m30 max
Step 55.0°C 0m30 max
Step 72.0°C 0m30 max
-----
Final Extn 72.0°C 5m
Final hold 4°C 10m
    
```

3-Step PCR with multiple 'Stages'

```

Heated Lid          105°C
Preheat Lid         on
Pause               off
-----
In denat 94.0°C    5min
Hot Start           off
----- Stage 1 -----
Number cycles       10
Step 94.0°C 0m30 max
Step 60.0°C 0m30 max
Step 72.0°C 0m30 max
----- Stage 2 -----
Number cycles       20
Step 94.0°C 0m30 max
Step 55.0°C 0m30 max
Step 72.0°C 0m30 max
-----
Final Extn 72.0°C 5m
Final hold 4°C 10m
    
```

PROGRAMMING

```

* * New Program * * [ ]
2 STEP TEMPLATE 501  [ ]
3 STEP TEMPLATE 502  [ ]
[  ↑ ↓ , ← to select ]
    
```

Create a new program by selecting ** New Program ** from the menu and press "Enter" ↵

```

UNNAMED #
Heated Lid          105°C
Preheat Lid         on
[  ↑ ↓ , ← to edit  ]
[  ▾ to save       ]
    
```

Give the program a name and edit any of the default fields as described above.

To add a stage and then edit it

```

In denat 94.0°C 5m00
Hot Start           off
-----
[  ← to insert stage ]
    
```

Use the "Down Arrow" ↓ key to scroll to the dotted line below 'Hot Start'; it will start to flash.

Press "Enter" ↵ to insert a stage.

```

----- Stage 1 -----
Number cycles          █
Step  94.0°C 00m30  max
[    ← edit no.cycles  ]

```

A new stage will be inserted which consists of a 'Number cycles' field and one step.

Edit the number of cycles by typing the number required then press "Enter" ↵

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C 00m30  max
[  '.' to insert step  ]
[    ↑ ↓, ← to edit   ]
[    ← to delete step ]

```

Use the "Down Arrow" ↓ key to move to the step.

Press "Enter" ↵ to edit the step as described above.

Press "Enter" ↵ again to accept.

To insert additional steps

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C 00m30  max
[  '.' to insert step  ]
[    ↑ ↓, ← to edit   ]

```

To insert a new step before the first step, use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to select 'Number cycles'.

Press the "Decimal Point" key to insert a new step.

To insert a new step after an existing step, use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to select the step.

Press the "Decimal Point" key to insert a new step.

To delete a step

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C 00m30  max
[  '.' to insert step  ]
[    ↑ ↓, ← to edit   ]
[    ← to delete step ]

```

Use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to select the step you wish to delete.

Press "Delete" ← to delete the step.

```

----- Stage 1 -----
Number cycles          10
Step  94.0°C 00m30  max
[    DEL STEP ?      ]
[    ← Yes or ▽ No   ]

```

The prompt line will flash between 'DEL STEP?' and 'YES or NO'.

Press "Enter" ↵ to delete the step.

or

Press "End" ▽ to return to editing the step.

To add a new stage before an existing one

```

----- Stage 1 -----
Number cycles           10
Step  94.0°C  00m30  max
[    ↵ to insert stage ]
[    ← to delete stage ]
    
```

Use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to highlight the stage number.

Press "Enter" ↵

A new stage will be added before the existing stage. The other stage numbers will be adjusted as appropriate. The new stage can now be edited as described above.

To add a new stage at the end

```

-----
Final Extn 72.0°C      5m00
Final hold           10.0°C
[    ↵ to insert stage ]
    
```

Use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to highlight the dotted line below the last stage.

Press "Enter" ↵

A new stage will be added with the next consecutive number. The new stage can now be edited as described above.

To delete a stage

```

----- Stage 1 -----
Number cycles           10
Step  94.0°C  00m30  max
[    ↵ to insert stage ]
[    ← to delete stage ]
    
```

Use the "Up Arrow" or "Down Arrow" ↑ ↓ keys to highlight the stage you wish to delete.

Press "Delete" ←

```

----- Stage 1 -----
Number cycles           10
Step  94.0°C  00m30  max
[    DEL STAGE?      ]
[    ↵ Yes or ▽ No  ]
    
```

The prompt line will flash between 'DEL STAGE?' and 'YES' or 'NO'.

Press "Enter" ↵ to delete the stage.

To insert a 'pause' into a program

```

Step  94.0°C  00m30  max
----- Stage 2 -----
Number cycles  █
Step  94.0°C  00m30  max

[  '.' to insert step  ]
[  ↑ ↓ , ↵ to edit    ]
    
```

To insert a pause after an existing stage, follow the steps for inserting a stage as above.

Highlight the 'Number cycles' line then press "Pause" | |

```

Step  94.0°C  00m30  max
----- Stage 2 -----
Number cycles  0
Step  94.0°C  00m30  max

[  CHANGE TO PAUSE    ]
[  ↵ Yes or ▽ No     ]
    
```

The prompt line will flash between 'CHANGE TO PAUSE?' and 'YES or NO'.

Press "Enter" ↵ to insert a pause.

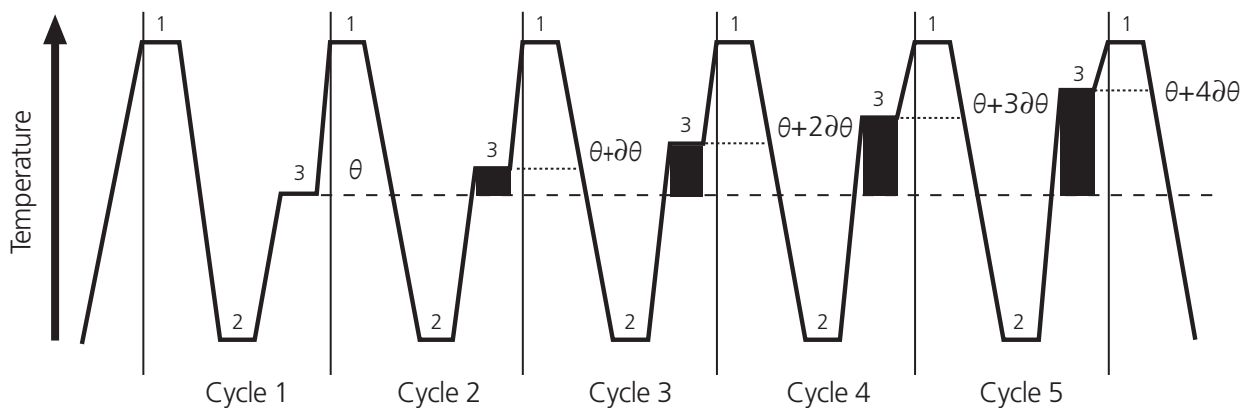
```

----- Stage 2 -----
Pause
-----
[  ↵ to insert step  ]
[  ← to delete step  ]
    
```

The stage number will then be highlighted with 'Pause' as the only line.

TO PROGRAM INCREMENT OR DECREMENT TEMPERATURE OR TIME


Under normal circumstances, the Hold Temperature of all steps is constant. However, it is possible to automatically increment or decrement the temperature of a specified step of a programme.

Temperature profile of incremented Hold Temperature at step 3 of each cycle

The Hold Temperature of the incremented/decremented step is the defined Hold Temperature plus the summation of the increments/decrements. The first cycle is never incremented/decremented, only subsequent cycles. If you select a decrement, the Hold Temperature is prevented from falling below 4°C. If you select an increment, the Hold Temperature is prevented from rising above 99°C.

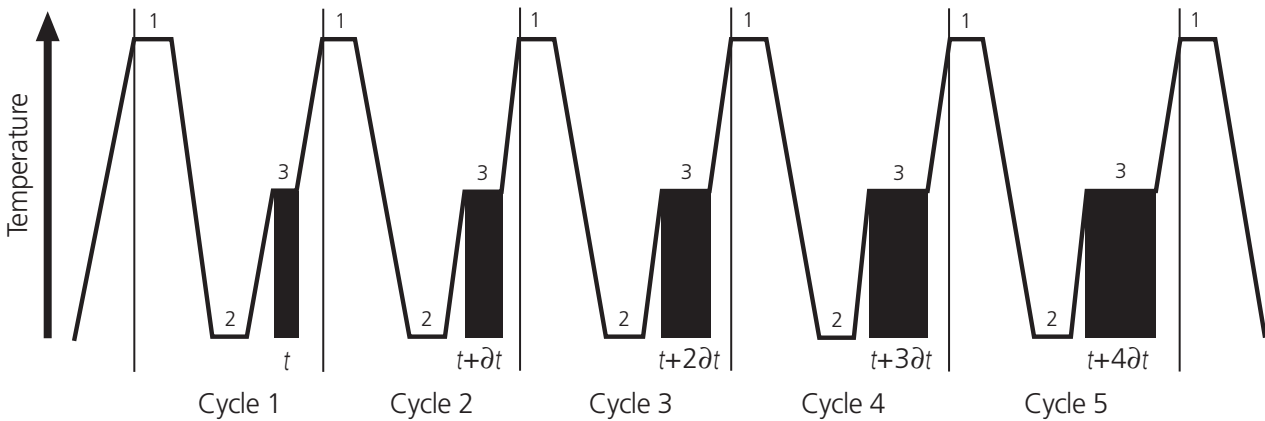Similarly, under normal circumstances, the Hold Time of all steps is constant. However, it is possible to automatically increment or decrement the duration of a specified step of a program.

Temperature profile of incremented Hold Time at step 3 of each cycle

The Hold Time of the incremented/decremented step is the defined Hold Time plus the summation of the increments/decrements. The first cycle is never incremented/decremented, only subsequent cycles. If you select a decrement, the Hold Time is prevented from falling below one second.

```

Number cycles          30
Step  94.0°C  00m30  max
Step  58.0°C  00m30  max
[  ↑ ↓ , ← to edit  ]
    
```

```

[  ← to delete step  ]
[  `.' to insert step ]
    
```

```

Number cycles          30
Step  94.0°C  00m30  max
Step  58.0°C  00m30  max
[  CHANGE STEP ?    ]
    
```

```

[  ← Yes or ▽ No  ]
    
```

Use the "Up Arrow" or "Down" Arrow" ↑ ↓ keys to move to the step where you want an increment or decrement temperature or time.

Press "Enter" ← to go into edit mode.

Press "Pause" | |

The prompt line will change to flash between 'CHANGE STEP?' and 'YES' or 'NO'.

Press "Enter" ←

The cursor will again move into the edit mode.

Now press the "Up Arrow" key ↑

```

Number cycles          30
Step  94.0°C  00m30  max
█ fst  58.0°C  00m30  max
[   ↑ ↓, ← to edit   ]
[       ▽ to save    ]
    
```

The step will change to 'fst', meaning "first", and the software will insert another line after it entitled 'lst', meaning "last".

Press "Enter" ↵ twice to go into edit mode.

```

Number cycles          30
Step  94.0°C  00m30  max
  fst  65█ °C  00m30  max
[   edit temperature  ]
    
```

Use the "Alpha/numeric" keys to change the values of the 'fst' line to the first temperatures and/or times you require.

Press "Enter" ↵ after each field you change.

You will exit 'Edit Mode' when you reach the end of the line.

```

Step  94.0°C  00m30  max
fst   65.0°C  00m30  max
lst  55.0°C  00m30
[   ↑ ↓, ← to edit   ]
[ ← to delete step   ]
[ `.' to insert step ]
    
```

Use the "Down Arrow" key ↓ to move onto the 'lst' line.

Press "Enter" ↵ to edit the step.

```

Step  94.0°C  00m30  max
fst   65.0°C  00m30  max
lst  55.0°C  1m30█
[   edit time        ]
[   . For hrs/mins   ]
    
```

Use the "Alpha/numeric" keys to change the values of the 'lst' line to the final temperatures and/or times you require.

Press "Enter" ↵ for each field you change.

You will exit 'Edit Mode' when you reach the end of the line.

When you run the program it will automatically increment or decrement in equal steps from the first time and/or temperature to the last over the number of cycles in the stage. For example: if the first temperature is 65°C and the last temperature is 55°C, over 30 cycles the temperature will decrease 0.33°C each cycle. Similarly, in the example shown, the time would increase from 30 seconds to 1 minute 30 seconds in 2 second increments.

To save the program

```

----- Stage 1 -----
Number cycles           10
Step  94.0°C   00m30   max
[                SAVE?                ]

[   ← Yes or ▽ No   ]

```

```

                GEORGE      (10)
Run Program
View Program
[   ↑ ↓ , ← to select   ]

```

A program can only be run once it has been saved. Press the "End" ▽ key.

The prompt line will flash between 'SAVE?' and 'YES or NO'.

Press "Enter" ↵ for 'Yes' to save the changes you have made.

If you press "End" ▽ for 'No' then all the changes you have made will be lost.

The program is now ready to run and will be allocated with a unique "Speed dial" number which can be used to rapidly locate the program from the main start-up screen.

Use the "Up Arrow" and "Down Arrow" ↑ ↓ keys to scroll through the list of options.

Stopping what you are doing

At any time during programming the TC-412, when you are satisfied with the program, you can finish and accept what you have done by first pressing the "End" ▽ key.

Then follow the instructions on screen. If you are not sure, wait for the instructions in the prompt line to scroll back to the first line that you see.

TO RUN A PROGRAM

```

Programs [0]
Information
Configuration
[ ↑ ↓ , ← to select ]
    
```

```

Programs [7]
Information
Configuration
[ ↑ ↓ , ← to select ]
    
```

```

* * New Program * * [ ]
2 STEP TEMPLATE 501 9
3 STEP TEMPLATE 502 9
[ ↑ ↓ , ← to select ]
    
```

```

DAVID 8 [D]
DONNA 7
[ ↑ ↓ , ← to select ]
    
```

```

DONNA [7]
Run Program
View Program
[ ↑ ↓ , ← to select ]
    
```

```

DONNA

[ RUN PROGRAM? ]

[ ← Yes or ▽ No ]
    
```

When the unit is switched on the display will show this screen.

If known, type in the 'Speed dial' number for the required program and press "Enter" ← This will take you directly to the required program.

If you do not know the 'Speed dial' number, press "Enter" ← to access the list of stored programs.

Next either:

1. Scroll down to the required program using the "Up Arrow" and "Down Arrow" ↑ ↓ keys.
2. Or type the first letter of the program name in the "New program" field; this will then list all the program names beginning with that letter. Use the "Up Arrow" and "Down Arrow" keys ↑ ↓ to select the required program.

Press "Enter" to accept ← For example 'Donna'.

To run the program, select 'Run program' and press "Enter" ←

A screen similar to this one will appear.

Press "Enter" ← to start the run.

```
Cycle 4/30          56m08
Sample  58.0°C
Hold @  58.0°C      00m17
[      ▽ to STOP   ]

[      || to PAUSE ]
[      ↵ for info  ]
```

During the run the LCD screen will report:

1. The cycle number
2. The time remaining (counting down in seconds)
3. The sample temperature (whether holding or ramping)
4. The current step of the run
5. A count down of the hold time.

An example is shown here.

Manual pause

```
Cycle 24/30        19m26
Sample  72.0°C
Hold @  72.0°C     00m10
[      PAUSE?     ]

[      ↵ Yes or ▽ No ]
```

To pause a program while it is running, press "Pause" || The run will continue but the prompt line will alternate 'PAUSE?' and 'YES or NO'

To confirm that you want the program to pause press "Enter" ↵

If you have pressed the "Pause" key by mistake and you want the program to continue, press "End" ▽

The unit will emit a beeping sound while paused.

To resume the run, press "Pause" || followed by "Enter" ↵ to continue.

```
Cycle 24/30        19m26
Sample  72.0°C
Paused
[      ▽ to STOP   ]

[      || to CONTINUE ]
[      ↵ for info    ]
```

Manual stop

```
Cycle 24/30        10m39
Sample  94.0°C
Hold @  94.0°C     00m10
[      STOP PROGRAM? ]

[      ↵ Yes or ▽ No ]
```

To stop a program while it is running, press "End" ▽

The run will continue but the prompt line will alternate 'STOP PROGRAM?' and 'YES or NO'

To confirm that you want to stop the program press "Enter" ↵

If you have pressed the "End" ▽ key by mistake and you want the program to continue, press "End" ▽ again.

When a program is stopped the unit will emit a beeping sound and the original program screen will appear.

Program Information

```

DONNA
Heated Lid           on
Status              ramping
[   ← to exit   ]
    
```

Pressing "Enter" ← while a program is running will give information about the current run including which program is actually running and the unit status e.g. holding or ramping.

An example is shown here.

INFORMATION

```

Programs             [0]
Information
Configuration
[   ↑ ↓, ← to select   ]
    
```

The **Information** field on the front screen contains all the unit-specific information such as serial number and software version.

Press "Enter" ← to access the unit information.

```

Block                96x0.2ml
Block s/n            139160-2
Unit s/n             131625-1
S/w version          33.01
H/w version          1
Comms address        1
Unit cycles          624
Block cycles         69
    
```

The complete list is displayed here.

Note that the "Comms address" must be unique for each unit if more than one cyclor is linked to a PC using the Gensoft control software.

CONFIGURATION

Warning! The Configuration software should only be accessed by designated personnel, as some parameters can be changed or may be lost when using it.

```

Programs             [0]
Information
Configuration
[   ↑ ↓, ← to select   ]
    
```

The **Configuration** field on the front screen allows a system administrator access to set up all the default settings for the programs, time and date settings and to change and remove passwords.

```

CONFIGURATION

Password            pad█
[   ← to accept   ]

[   ▽ to exit   ]
    
```

Access to the 'Configuration' set up is password protected.

To enter configuration, use the alpha keys to type in 'pad'

Press "Enter" ←

```

Program defaults
Settings
Passwords
Controller reset
    
```

There are a number of options which can be changed; these are listed here.

Use the "Up Arrow" and "Down Arrow" ↑ ↓ keys to select then press "Enter" ↵ to access any of the options. Each option is discussed in detail below.

Program defaults

```

Program defaults
Settings
Passwords
Controller reset
    
```

Highlight 'Program defaults' then press "Enter" ↵

```

Heated Lid           105°C
Preheat Lid           off
Pause                 off
-----
In denat              94°C      5m00
Hot Start              off
-----
Final Extn            72.0°C    5m00
Final hold             10.0°C
-----
Lid type               grey
[   ↑ ↓ , ↵ to edit   ]
    
```

The complete list of program defaults is given here. These are the values that all new programs will have when you start to create a new program.

To edit any line use the "Up Arrow" or "Down Arrow" keys ↑ ↓ to select the line you wish to change.

Press "Enter" ↵ then change the value using the numerical keys. Press "Enter" ↵ to accept the new value.

To edit a field with on/off settings, first select the field and press "Enter" ↵

Use the "Up Arrow" and "Down Arrow" ↑ ↓ keys to toggle between the settings. Press "Enter" ↵ to accept.

When you have set all the defaults, press the "Enter" ↵ key to go back to the Configuration menu.

Note: altering the default values does not change existing programs.

Settings

```

Program defaults
Settings
Passwords
Controller reset
    
```

Highlight 'Settings' then press "Enter" ↵

```

Set date      xx / xx / xx
Set time      xx : xx
Date format   dmy
Auto restart  on
Beeper        on
Coms address  1
Coms speed    4800
[  ↑ ↓ , ← to select  ]
    
```

The complete list of settings is given here.

Auto restart "on" will allow the unit to restart from where it left off should there be a power failure and the power returns.

The "Comms address" refers to the location of a unit in a daisy chain and must be unique if there are several units linked to each other.

The optimum "Coms speed" is set at 4800. When using the Gensoft control software, ensure all cyclers are set to the same value.

Edit any line as described for 'Program defaults' above. Press "Enter" ← to accept the new value.

When you have finished, press the "Enter" ← key to go back to the 'Configuration' menu.

Passwords

```

Program defaults
Settings
Passwords
Controller reset
    
```

The 'Passwords' function allows the administrator to change the 'Admin password' (which allows access to Configuration) and also to delete any passwords associated with programs.

Highlight 'Passwords' then press "Enter" ←

```

* * Admin passwd * *
DONNA
GEORGE
[  ↑ ↓ , ← to select  ]
    
```

The screen will appear as opposite.

We suggest that you change the 'Admin password' at this stage in order to protect the configuration settings.

```

      ADMIN PASSWD
Password  █
Confirm
[  ← to accept  ]
    
```

Use the "Up Arrow" ↑ key to highlight 'Admin passwd

'Press "Enter" ← and type in the new password. Repeat to confirm and accept.

The screen will go back to the 'Configuration' menu.

```

* * Admin passwd * *
DONNA
GEORGE
[  DEL PASSWORD?  ]
[  ← Yes or ▽ No  ]
    
```

To delete a password associated with a program, use the "Up Arrow" and "Down Arrow" ↑ ↓ keys to highlight the program.

Press "Enter" ←

Press "Enter" ← again to delete the password.

Controller reset

```

Program defaults
Settings
Passwords
Controller reset
    
```

The 'Controller reset' will delete all saved programs from the unit.

Highlight 'Controller reset' then press "Enter" ↵

```

          CONTROLLER RESET
This will delete
all programs
[          SURE ?          ]
    
```

To delete all the programs, press "Enter" ↵

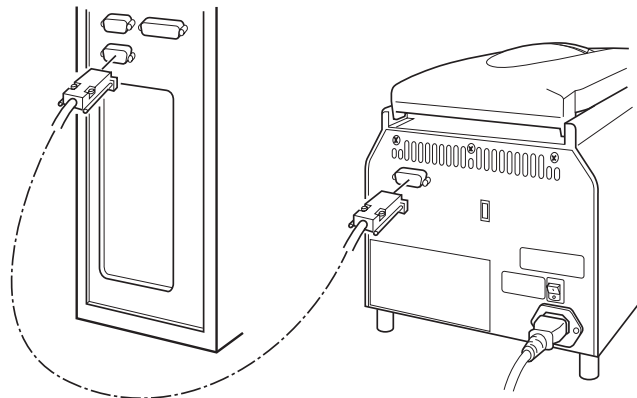
```

[          ↵ Yes or ▽ No          ]
    
```

RUNNING THE TC-412 FROM A PC

The TC-412 can be run from a PC using Techne's Gensoft software. This is available to download from the Techne web-site at www.techne.com. You will need 'Winzip' (which is available free of charge on www.winzip.com) to unzip the files. Instructions for using Gensoft are included in a separate manual.

CONNECTING ONE UNIT TO THE COMPUTER USING AN RS232 CABLE

If you are driving one TC-412 unit you will need part number FGEN232 which includes a cable, Gensoft on disk and an instruction sheet.


CONNECTING SEVERAL UNITS TO THE COMPUTER USING THE DAISY CHAIN CABLES

If you are driving two or more units then you will need part number FGEN485D, a 230V UK power pack set, or FGEN485E, a 230V European power pack set, or FGEN485P, a 120/100V USA power pack set. A power pack set includes: the connection to a computer (which also converts to RS485), a power pack, the necessary cables, a termination box and the software on disk with an instruction sheet. You will also need one or more of the following:

-
- FAGENTWO** A two unit set which includes two connectors for the cyclers and a cable to interconnect them.
 - FGENFIVE** A five unit set which includes five connectors for the cyclers and four cables to interconnect them.
 - FAGENTEN** A ten unit set which includes ten connectors for the cyclers and nine cables to interconnect them.
 - FGENONE** A one unit set for extending the daisy chain further which includes one connector for the cyclers and one cable for connecting it to the daisy chain.
-

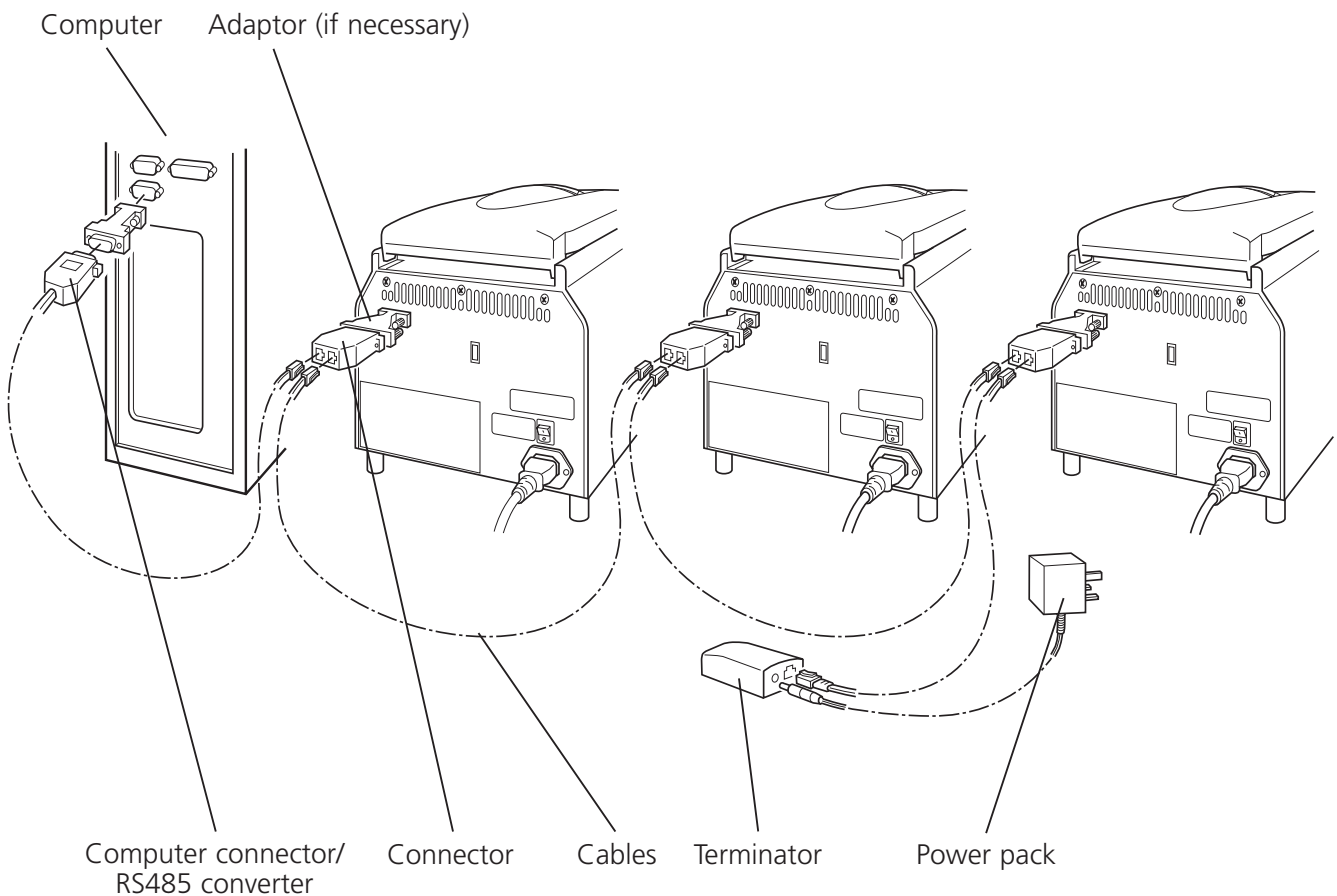Notice from the diagram that the cable connections to the cyclers are alternately bottom to top on the connector.

UPGRADING THE SOFTWARE IN YOUR TC-412

Connect your TC-412 to your computer using an RS232 lead. Log onto the Techne web-site at www.techne.com. Click on "Thermal Cycler Software Update" at the top of the page and you will find information on how to upgrade your software. Click on 'Update Software' and follow the instructions on the page to install the programs.

[SEARCH]

USER MAINTENANCE

CLEANING YOUR TC-412

Before cleaning your unit, disconnect from the power supply and allow to cool to below 50°C. The heating/cooling block, including wells and flat surfaces, should be cleaned regularly to ensure optimum heat transfer to the samples. Always clean the block if there has been a spillage. Use a cloth or cotton buds dipped in a fresh, 50:50 water/isopropanol solution and make sure that no deposits are left in the wells.

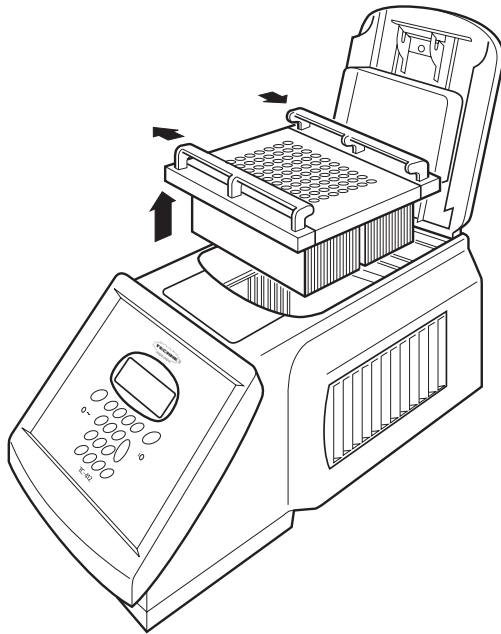
In the case of radioactive spillages the heater block can be removed from the unit for more detailed cleaning. Techne recommends that you use a proprietary cleaning agent and follow the manufacturer's instructions. The heating/cooling block is made of aluminium, therefore an agent such as Neutracon (from Decon Labs Ltd.), suitable for nonferrous metals should be used. However, remember that other parts of the unit are made of ferrous materials and may be damaged by spillage onto them.

The case of the TC-412 may be cleaned with a cloth dipped in water or ethanol (methanol or formaldehyde can also be used). No part of the case or cover should be immersed in the solvents. Do not use aggressive solvents such as acetone or abrasive cleaners.

Before using any cleaning or decontamination method except those recommended here, the responsible person should check with Techne that the proposed method will not damage the equipment.

BLOCK REMOVAL

Before removing the block, disconnect from the power supply and allow it to cool to below 50°C.

- Slide the quick release handles in the direction of the arrows indicated and lift the block assembly out of the unit.
- Never lift or carry the block by one quick release handle, always use both handles or support the block from underneath.
- When re-fitting the block, or fitting a new one, slide the quick release handles in the opposite direction to lock the assembly into the unit.
- Open or closed padlock symbols are printed on the quick release handles for easy reference.

FREQUENTLY ASKED QUESTIONS

Q1 What is the 'Pause' function at the start of the program used for?

A1 Some users prefer to preheat the heated lid before placing the samples into the unit. The pause feature is used to stop the unit after the 4 minute heated lid preheat step. It will also sound an audible alarm indicating that the machine is ready for the sample tubes or plate to be added. Pressing the "Pause" key followed by the "Enter" key will commence the remaining program.

Q2 Why do you have to press some keys twice to make the unit do what you want it to do?

A2 With many of the program features there is a fail-safe function which asks if you are sure you want to perform a particular action. You are required to press one key to select a function followed by a second to accept it.

Q3 What is 'Hot Start'?

A3 A "hot start" is used for reducing non-specific product formation during the first cycle of amplification. The 'Hot Start' programming step is used to pause the instrument at a specific temperature, typically around 70°C, after the initial template denaturation. The reason is to allow the manual addition of unmodified Taq DNA polymerase which may lose activity if added during the initial 5min denaturation. Heat-activated Taq or Hot Start enzymes do not require this step.

Q4 What is the incremental function?

A4 Incremental timing and temperature are used to increase or decrease either the time or temperature incrementally over the number of cycles in a stage.

Incrementation of extension time is used with 'Long PCR' which is when large template fragments are to be amplified (e.g. 27kb lambda DNA, 40kb genomic DNA).

Decremental temperature is used for protocols such as 'Touchdown PCR' where one starts with a high annealing temperature in the first cycle and gradually decreases the temperature over the number of cycles in the stage. This ensures that only the specific product is amplified.

Q5 What is the appropriate pressure of the heated lid?

A5 The lid will need to be adjusted for the height of each different type of consumable used in the unit. Firstly, with the lid in the down position, turn the orange rotating knob anticlockwise until it will not go any further. Next with the tubes or plate in the unit, lock the lid and turn the knob clockwise until pressure on the tubes or plate is just felt; then rotate another quarter turn. Do not over-tighten.

If you require further technical or applicational assistance please contact Techne at:
technehelp@barloworld-scientific.com.

For servicing information please contact: **service@barloworld-scientific.com.**

We are continually striving to improve our thermal cyclers and software. If you have any comments and suggestions on how we can do things better please send them to us at:
techne@barloworld-scientific.com

ADDITIONAL INFORMATION

Brief fault finding notes and a list of replacement parts are given in this section.

Note that this equipment should only be dismantled by properly trained personnel. Removing the outer cover exposes potentially lethal mains voltages. There are no user serviceable parts within this equipment.

FAULT FINDING

Should you have any problems with your TC-412 which cannot be easily remedied, you should contact your supplier and return the unit if necessary. Please include details of the fault observed and remember to return the unit in its original packing. Techne accepts no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier, giving the full serial number of the unit and software version number (shown when the unit is first switched on).

FUSES

If neither the power light nor display on the front panel is lit, one of the two fuses may have blown. Check that there is no external cause, such as a faulty plug or lead. Check both fuses and replace the faulty fuse with a new one of the correct value (fuse values are given on the label next to the power inlet). Note that fuses should only be replaced by a qualified electrician.

The holder for the two fuses is built into the mains input socket. First remove the power cable and then gently prise the fuse drawer open with a flat-bladed screwdriver or similar tool.

Each fuse can be removed by using the screwdriver as a lever.

Exchange the faulty fuse in the fuse holder for a working fuse of the correct value. Finally, replace the fuse drawer in the fuse compartment and push the drawer shut.

Fuses which blow repeatedly indicate a serious fault and you should return the unit to your supplier for repair.

THE HEATED LID OVER-TEMPERATURE CUT-OUT

The heated lid is fitted with an independent circuit to protect it from overheating. In the unlikely event of an over-temperature problem with the lid, the unit is fitted with a thermal fuse which will remove power to the heater plate should the maximum temperature be exceeded.

INSULATION TESTING

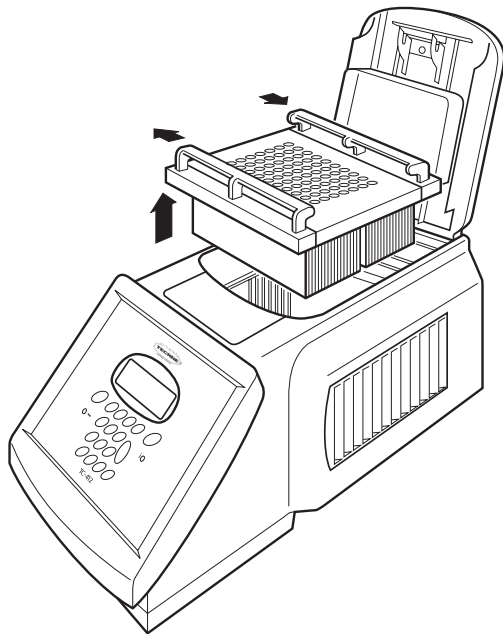
This equipment is fitted with RFI suppression circuitry. Any check of the electrical insulation by means of high voltage dielectric testing (for example as in BS EN 61010-1) must be carried out using only a DC voltage.

This unit contains semiconductor components which may be damaged by electric field effects.

INTERCHANGEABLE BLOCKS

The block can be removed and replaced by another of the same sort or of a different sort. The software will register the change and set up the unit for the new block.

Before removing the block, disconnect from the power supply and allow it to cool to below 50°C.

- Slide the quick release handles in the direction of the arrows indicated and lift the block assembly out of the unit.
- Never lift or carry the block by one quick release handle, always use both handles or support the block from underneath.
- When re-fitting the block, or fitting a new one, slide the quick release handles in the opposite direction to lock the assembly into the unit.
- Open or closed padlock symbols are printed on the quick release handles for easy reference.
- Remember to adjust the options in the Options menu for the new block.

If you have fitted a flat block for slide applications, you must set the heated lid to OFF.

ACCESSORIES

The following accessories can be obtained from Techne or your Techne distributor:

Item No	Description	Quantity
FTC41B5D	60 x 0.5ml micro tubes	1
FTC41BHD	96 x 0.2ml micro tubes or 96 well plate	1
FTC41BFD	96 x 0.2ml micro tubes or 96 well fully skirted plate	1
FTC4B384	384 well block	1
FTC41BID	Flat plate (<i>in situ</i>) block	1
FTUB02TW	0.2ml micro tubes	Pack of 1000
FTUB05TW	0.5ml micro tubes thin wall	Pack of 1000
FMIP384	384 well plate	Pack of 50

REPLACEMENT PARTS

The following replacement parts can be obtained from Techne or your Techne distributor:

Item No	Description	Quantity required
FCABRTUK	Mains cable and plug, 230V UK	1
FCABRTEU	Mains cable and Schuko plug, 230V Europe	1
FCABRTUS	Mains cable and 3-pin plug, 100V/120V US type	1
6500342	Fuse T6.3A (230V units)	2
6008445	Fuse T8A (120/100V units)	2
6500324	Fuse T10A (100V units)	2

**Barloworld
Scientific**

Barloworld Scientific Ltd

Beacon Road Stone
Staffordshire ST15 0SA
United Kingdom
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 813748
e-mail: techne@barloworld-scientific.com
www.barloworld-scientific.com

Barloworld Scientific US Ltd

Techne
3 Terri Lane Suite 10
Burlington N.J. 08016, USA
Tel: 800-225-9243
Fax: 609-589-2571
e-mail: sales@techneusa.com
www.techneusa.com

Barloworld Scientific France SAS

ZI du Rocher Vert - BP 79
77793 Nemours Cedex
France
Tel: +33 1 64 45 13 13
Fax: +33 1 64 45 13 00
e-mail: bsf@barloworld-scientific.fr
www.barloworld-scientific.com

Barloworld Scientific Italia Srl

Via Alcide de Gasperi 56
20077 Riozzo di Cerro al Lambro
Milano Italia
Tel: +39 (0)2 98230679
Fax: +39 (0)2 98230211
e-mail: marketing@barloworld-scientific.it
www.barloworld-scientific.it